

Instruction Manual



Biosciences

LI-840A CO₂/H₂O Analyzer

Instruction Manual



Federal Communications Commission

Radio Frequency Interference Statement

WARNING: This equipment generates, uses, and can radiate radio frequency energy and if not installed in accordance with the instruction manual, may cause interference to radio communications. It has been tested and found to comply with the limits for a Class A computing device pursuant to Subpart J of Part 15 of FCC rules, which are designed to provide a reasonable protection against such interference when operated in a commercial environment. Operation of this equipment in a residential area is likely to cause interference in which case the user, at his own expense, will be required to take whatever measures may be required to correct the interference.

		LI-8	340A			
			Hazardous Su	ibstances or Ele	ements	
Component Name	Lead (Pb)	Mercury (Hg)	Cadmium (Cd)	Chromium VI Compounds (Cr ⁶⁺)	Polybrominated Biphenyls (PBB)	Polybrominate Diphenyl Ethe (PBDE)
Optical Bench Assembly	X	0	0	0	0	0
Source Block Assembly	х	0	0	0	0	0
Detector Block Assembly	Х	0	0	0	0	0
Ribbon Cable Assembly	0	0	0	0	0	0
Serial Cable Assembly	х	0	0	0	0	0
LED Circuit Board Assembly	0	0	0	0	0	0
RS-232 Cable Kit	0	0	0	0	0	0
Main Board V2 Circuit Board Assembly	Х	0	0	0	0	0
Red Cap Plug	0	0	0	0	0	0
X: this component does contain this haza	ardous subs	tance above th	e maximum cono	centration value	is in nomogeneou	is materials
specified in the SJ/T 11363-2006 Industr	y Standard	(Company can	explain the tech		or the "X")	
specified in the SJ/T 11363-2006 Industr	y Standard	(Company can	explain the tech		or the "X")	
specified in the SJ/T 11363-2006 Industr	y Standard	(Company can	explain the tech		or the "X")	
specified in the SJ/T 11363-2006 Industr			explain the tech	nical reasons fo	or the "X")	
specified in the SJ/T 11363-2006 Industr			H ₂ O 分析仪	nical reasons fo	or the "X")	
specified in the SJ/T 11363-2006 Industr 部件名称			H ₂ O 分析仪	nical reasons fo	or the "X") 多溴联苯 (PBB)	多溴二苯酮 (PBDE)
	LI-8	340A CO ₂ /	/H ₂ O 分析仪 _{有毒有}	nical reasons fo		
部件名称	LI-{ 铅 (Pb)	340A CO ₂ / 汞 (Hg)	/H ₂ O 分析仪 _{有毒有} 镉 (Cd)	nical reasons fo 害物质或元素 六价铬 (Cr ⁶⁺)	多溴联苯 (PBB)	(PBDE)
· 部件名称 光路系统组件	LI-8 倍 (Pb) ×	340A CO ₂ / 汞 (Hg) o	H ₂ O 分析仪 _{有毒有} 镉 (Cd) o	nical reasons fo 害物质或元素 六价铬 (Cr ⁶⁺) O	多溴联苯 (PBB) O	(PBDE) O
部件名称 光路系统组件 红外光源组件	LI-{ 铅 (Pb) X X	340A CO₂/ 汞 (Hg) 0 0	H ₂ O 分析仪 有毒有 锅 (Cd) 0	nical reasons fo 害物质或元素 六价铬 (Cr ⁶⁺) O	多溴联苯 (PBB) 0 0	(PBDE) O O
部件名称 光路系统组件 红外光源组件 红外光波调器组件	LI-{ 铅 (Pb) X X X	340A CO₂/ 汞 (Hg) ○ ○	H ₂ O 分析仪 有毒有 镉 (Cd) 0 0	nical reasons fc 害物质或元素 六价铬 (Cr ⁶⁺) 〇 〇 〇	多溴联苯 (PBB) 0 0 0	(PBDE) 0 0 0
部件名称 光路系统组件 红外无波组件 红外无论调器组件 带状也缆出件	LI-8 铅 (Pb) X X X X O	340A CO₂/	H ₂ O 分析仪 有毒看 锅 (Cd) 0 0 0	nical reasons fc 言物质或元素 六价铬 (Cr ⁶⁺) 〇 〇 〇	多溴联苯 (PBB) の の の	(PBDE) 0 0 0 0
部件名称 光路系线组件 红外无路组件 红外无路器组件 带状电缆组件 带状电缆组件	LI-{ 铅 (Pb) X X X X 0 X 0 0 0	340A CO₂/	H ₂ O 分析仪 有毒	nical reasons fc 害物原或元素 六价格 (Cr ⁶⁺) 0 0 0 0 0 0 0 0 0 0 0	多湊联苯 (PBB) 0 0 0 0 0	(PBDE) 0 0 0 0 0 0 0 0 0
部件名称 部件名称 近外光波调件 近外光波调告 近外光波调告 近外光波调告 重大电缆调件 带打电缆组件 每行电缆组件 每行电缆组件 每行电缆组件 专行影响。 2000 1	LI-{ 铅 (Pb) X X X X O X O	340A CO₂/ 〒 ○	H ₂ O 分析仪 有毒有	nical reasons fo 書物质或元素 六价铬 (Cr ⁶⁺) の の の の の の	多決联苯 (PBB) 0 0 0 0 0 0 0	(PBDE) 0 0 0 0 0 0
部件名称 光路系统组件 红外光波照件 证外光检测器组件 带状电缆组件 串行电缆组件 串行电缆组件 串行电路板组件 串行 RS-232 电缆	LI-{ 铅 (Pb) X X X X 0 X 0 0 0	340A CO₂/ (Hg) 0 0 0 0 0 0 0 0 0 0	H ₂ O 分析仪 有毒有 镉 (Cd) 0 0 0 0 0 0	nical reasons fc 害物原或元素 六价格 (Cr ⁶⁺) 0 0 0 0 0 0 0 0 0 0 0	多溴联苯 (PBB) 0 0 0 0 0 0 0 0	(PBDE) 0 0 0 0 0 0 0 0
部件名称 部件名称 近外光波调件 近外光波调告 近外光波调告 近外光波调告 重大电缆调件 带打电缆组件 每行电缆组件 每行电缆组件 每行电缆组件 专行影响。 2000 1	LI-{ 铅(Pb) X X X X O X O O X	★ (Hg) 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0 0	H ₂ O 分析仪 石毒石 幅 (Cd) 0 0 0 0 0 0 0 0 0 0 0 0 0	nical reasons fo 言物质或元素 六价铬 (Cr ⁶⁺) 〇 〇 〇 〇 〇 〇 〇 〇 〇 〇	多凍联苯 (PBB) 0 0 0 0 0 0 0 0 0	(PBDE) 0 0 0 0 0 0 0 0 0 0
部件名称 部件名称 近外光波调件 近外光波调告 近外光波调告 近外光波调告 重大电缆调件 带打电缆组件 每行电缆组件 每行电缆组件 每行电缆组件 专行影响。 2000 1	LI-{ 铅 (Pb) X X X X O X O X O X O O X	₩ (Hg) 0	H ₂ O 分析仪 有毒有 镉 (Cd) 0 0 0 0 0 0 0 0 0 0	ical reasons for 害物质或元素 六价铬 (Cr ⁵⁺) 0 0 0 0 0 0 0 0 0 0 0 0 0		(PBDE) 0 0 0 0 0 0 0 0 0 0

Doc. #53-13776-A

May 21, 2013 Page 1 of 1

LI-COR® Biosciences		LI-COR, inc. Environmental 46/7 Superior Street P.O. Box 4425 Lincoln, Netwaska 68504 USA Phone: 402-467-3576 FAX: 4024-6723174-5756 (U.S. & Cunada) Franal: envsales@licor.com
Ι	Declaration of Con	formity
Manufacturer's Name:	LI-COR Inc.	
Manufacturer's Address:	4647 Superior Street Lincoln, Nebraska USA 68504	L .
This declaration of conformit	y is issued under the sole resp	onsibility of the manufacturer.
Product Description:	The LI-840A is a CO2 and H2	O Gas Analyzer.
Product Name: Model Number(s): Product Options:	Gas Analyzer LI-840A, LI-840-01, LI-840A None	-02
	ne requirements of the EMC D rries the CE marking accordin	irective 2004/108/EC, and the RoHS gly.
IEC 61000-4-2 : 2008 IEC 61000-4-3 : 2006/ IEC 61000-4-4 : 2012 IEC 61000-4-5 : 2005 IEC 61000-4-6 : 2008	iated Emissions, Class A ducted Emissions, Class A uipment for measurement, contr ESD, 4KV/8KV Contact/Air A1:2007/A2:2010 Radiated RF EFT/Burst Surge Conducted RF Immunity H-Field Immunity, 3A/m	
FCC 47 CFR Part 15.109 Radia FCC 47 CFR Part 15.107 Cond		
EMC performance can be affect	standards listed above is claimed ted by the complete installation	I for the non-OEM version of the LI-840A. therefore if using OEM versions (LI-840- re-qualify the complete system to the EMC

Directive.

May 21, 2013 Page 1 of 1

Document #53-13759-A

Dala

John Rada Engineering Manager

NOTICE

The information contained in this document is subject to change without notice. LI-COR MAKES NO WARRANTY OF ANY KIND WITH REGARD TO THIS MATERIAL, INCLUDING, BUT NOT LIMITED TO THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. LI-COR shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

This document contains proprietary information which is protected by copyright. All rights are reserved. No part of this document may be photocopied, reproduced, or translated to another language without prior written consent of LI-COR, Inc.

© Copyright 2010-2016, LI-COR, Inc. Publication Number 984-10690

Printing History

April 2010 January 2011 Updated May 2013; May 2015; June 2016

New editions of this manual will incorporate all material since the previous editions. Update packages may be used between editions which contain replacement and additional pages to be merged into the manual by the user.

The manual printing date indicates its current edition. The printing date changes when a new edition is printed. (Minor corrections and updates which are incorporated at reprint do not cause the date to change).

LI-COR, Inc. • 4421 Superior Street • Lincoln, Nebraska 68504 Phone: 402-467-3576 • FAX: 402-467-2819 Toll-free: 1-800-447-3576 (U.S. & Canada) e-mail: envsales@licor.com • www.licor.com

Table of Contents

Section 1. General Description

What's What	1-1
Features	1-3
Precautions	1-3
Getting Started Tutorial	1-4

Section 2. Power On

Power On	2-1
Alarms	2-1

Section 3. Operation

Installing the PC Communication Software on Your Computer	3-1
Setting the Communication Parameters	3-2
Cabling	3-2
RS-232 Output	3-2

Initial Setup	3-3
Using the Toolbar	3-7
Settings Window - Setting Operational Parameters	3-8
Options	
Enable Heater	3-10
Pressure Compensation	
Filter	3-10
DAC Output	
Converting Voltage Output to Cell Temperature	
Converting Voltage Output to Cell Pressure	3-13
Converting Current Output to ppm CO2 or mmol/mol H2O	3-14
CO2 Span Range	3-15
Alarms	
Using the Terminal Strip	
Charting Window	3-18
Series 1	
Series 2	
X-Axis Max	
Diagnostic Window	
Logging Data	
Start Logging	3-21
Stop Logging	
Pause Logging	

Options	3-22
Calibration Window - Setting the Zero and Span	
Two Point Span (Span 2 Tab)	
Considerations for performing a secondary span:	
Coefficients Page	
Advanced Page	3-32
Connection the LI-840A to the LI-1400 Datalogger	

Section 4. Theory of Operation

System Overview	4-1
Optical Bench System	4-2

Section 5. Maintenance

Cleaning the Optical Bench	5-1
Changing the Fuse	5-5
Note on Infrared Source Replacement	

Appendix A. Specifications

Appendix B. Equation Summary

General Form	. B-1
H2O Equations	
CO2 Equations	. B-3
Zeroing H2O	
Zeroing CO2	B-7
Spanning H2O	B-7
Secondary Span H2O	B-8
Spanning CO2	. B-9
Secondary Span CO2	B-10
Symbol Summary	B-10

Appendix C. Pin Assignments

DB-9 Connector

Appendix D. Suppliers

Appendix E. Configuration Grammar

Introduction	. E-1
LI-840A Communications and XML	
Connecting and Configuring Data Output	. E-3
Sending Data to the LI-840A	
Reading Data From the LI-840A	E-5
Polling the Current State of the LI-840A	E-5
Calibration	. E-6
To Zero the LI-840A	E-7
To Span the LI-840A	. E-8
To Perform a Two Point Span	. E-8
LI-840A XML Grammar and Element Description	E-10
Data Types in the XML Grammar	E-10
XML Grammar	

Warranty

General Description

What's What

This procedure should be followed if you have just taken delivery of your LI-840A. Check the packing list to verify that you have received everything that was ordered and that you have also received the following items:

- **RS-232 Cable** (Part #392-04993), used to connect the LI-840A to your computer. This cable is a null modem (cross-over) type cable; a straight through type cable will not work with the LI-840A. Appendix C shows the pin assignments for the DB-9 connector on the LI-840A front panel. An RS-232-to-USB converter is provided as well (Part # 6400-27)
- Data Communications Software This Windows® 7, 8/8.1, and 10 compatible software is used for communication between the LI-840A and a computer. The software is provided on CD. A complete description of this software can be found in "*Operation*" on page 3-1.

• Standard Spare Parts Kit (9980-021) - This kit contains replacement parts for the LI-840A. As you become familiar with the analyzer you will learn which items to keep close at hand and which items can be stored away. The spare parts kit contains these items:

Description	Qty.	LI-COR Part No.
Polyurethane Mounting Feet	4	234-02268
Mounting Screws	4	149-02610
Terminal Block	1	331-05273
Bev-a-line Tubing	12'	222-01824
Gelman Air Filter	2	9967-008
Quick Connect Unions	2	300-03123
Rapid Tube Fitting Nut	2	9861-036
2A Fuse	2	439-03952
Cleaning Kit, includes:	1	9980-013
5" Cleaning Swabs	5	610-05315
O-Rings	4	192-00226
Pre-formed Optical Bench Tubing	4	6580-010

Features

The LI-840A is an economical, high performance, non-dispersive infrared gas analyzer designed to be used for a wide variety of applications. Some of the LI-840A's important features include:

- A 14 cm (5.5") optical bench is used to obtain a CO₂ measurement range of 0-20,000 ppm (parts per million, or µmol/mol) and an H₂O measurement range of 0-80 mmol/mol.
- The optical bench is fully serviceable by the user; cleaning the optical bench is a simple process, and does not affect the instrument's factory calibration.
- Simple Windows® software provides for easy user calibration and data output.

Precautions

- The optical bench is maintained at a constant temperature of about 50 °C. Avoid direct exposure to sunlight or extremely high temperatures that may elevate the temperature inside the LI-840A case.
- The optical source is sensitive to vibration, and can be damaged by strong vibrations or jarring. Do not drop the LI-840A or expose it to severe mechanical shock.

- Do not use abrasive cleansers when cleaning the optical bench, as damage to the gold plating may result. Instructions for cleaning the optical bench can be found in Section 5.
- Always filter air entering the LI-840A. Two Gelman filter assemblies (p/n 9967-008) are included in the spares kit.

Getting Started Tutorial

The following section briefly covers the basic steps you might follow to set up the LI-840A to collect and record data. Many of these steps are described in greater detail throughout this manual.

Install the PC Communications Software.

This software is used to set the zero and span of the LI-840A and set up data communication and data transfer parameters for viewing and logging data. Installation instructions can be found in *"Installing the PC Communication Software on Your Computer" on page 3-1.*

2

Determine your jumper settings for high and low alarm output, if required.

There are 2 jumpers inside the LI-840A case (on the main circuit board), which toggle the high and low alarm output between 0-5V (TTL levels) and an open drain condition.


The default jumper positions are set for 0-5V output for both high and low alarms. This alarm configuration is suitable for logic devices such as dataloggers. To use the high and/or low alarms to operate a relay switch, the jumpers should be reconfigured for open drain output. See "Alarms" on page 2-1, for a description of how to change the alarm jumper settings.

If the default jumper positions are suitable for your application, or if you do not intend to use the alarms, proceed to Step 3.

3

Connect a power source to the LI-840A

The LI-840A requires an input voltage of 12-30 VDC. The power supply must be able to source a maximum current drain of 1.2A (at 12 VDC). After the instrument has warmed up it will draw about 0.3A (at 12 VDC) with the heaters on.

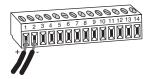

Bare wire leads are connected to the terminal strip at positions 1 and 2. Connect the positive lead (usually red) to position #1 (12-30VDC), and the negative lead to position #2 (GND).

NOTE: Pull straight out on the face of the terminal strip to remove it; this can make it easier to connect the wires.

4

Connect the RS-232 cable.

Connect one end of the serial cable to the 9-pin Serial I/O port on the LI-840A front panel, and the other end to a serial port on your computer. If you want to interface to a device with a 25-pin serial port, a 9-pin to 25-pin adapter must be used; a gender changer may be required in some cases. Tighten the two screws on the ends of the serial cable.

5


Connect other output device(s) if desired.

Output devices such as dataloggers and relay switches for high and low alarms are connected via the terminal strip. A complete description of the terminals can be found in Section 3, Using the Terminal Strip.

6

Connect an air pump and external power source.

The LI-840A sample cell must be continually flushed while making measurements, which requires an external air pump that supplies a flow rate of between 0.25 and 1.0 liters per minute. LI-COR offers a Diaphragm Pump under part number 286-04198 that is suitable for use with the LI-840A. One pump plumbing option is shown below.

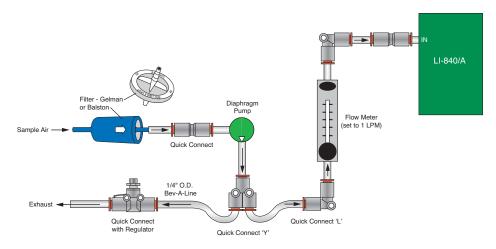




NOTE: More information about plumbing options and parts available from LI-COR can be found in Application Note #IRG4-105, entitled 'Air Pumps and Accessories for LI-COR Gas Analyzers - Suggested Parts and Vendors', which can be downloaded from www.licor.com/env/support.

Connect the input air stream and external filter.

Remove the nut from the **FLOW IN** fitting and conect a short (12" or so) piece of Bev-a-line tubing (from the spares kit). Tighten the nut. Place a Geman filter (also from the spares kit) in the sample airstream before it enters the **FLOW IN** port on the LI-840A. A user-supplied pump is required to force air through the LI-840A.

IMPORTANT: Always install the external air filter before operating the LI-840A. Insert filter into the input airstream before it enters the LI-840A. Failure to do this will lead to contamination of the optical path.

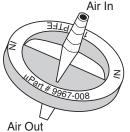

When using the LI-840A for applications where sample air is particularly dirty, you may consider stacking two filters in series. For maintenance, replace the filter furthest upstream from the analyzer with the filter closest to the analyzer and replace the filter closest to the analyzer with a new filter.

Install the new filter(s) with the blue lettering facing away from the FLOW IN port.

The air inlet and outlet ports should be covered with the dust caps provided or attached to an air line when the instrument is not in use. This will prevent dust from entering the instrument downstream from the filters where it can enter the optical path.

Replacement Gelman Acro 50 filters can be obtained from LI-COR (part #9967-008).

8

Start the PC Communications Software.

Click on the program icon to start the application. The Main window appears. Select **Connect** from the File menu. Choose the serial port to which the LI-840A is connected, and the data output interval. Click the **Connect** button. If the analyzer is connected properly, data will appear in the main window.

9

Configure the analyzer.

Select **Settings** from the View menu. The Settings window appears. Choose a value for signal averaging (the Filter field), between 0 and 20 seconds. Set values for high and low alarms and corresponding dead bands, if desired. See Section 2 for a complete description of the alarms. If you are using an output device to collect data, choose the value to be output at DAC 1 and/or DAC 2. Make sure the Enable Heater and Pressure Compensation buttons are enabled (checked), and click **OK**.

10 Set the analyzer zero. NOTE: The LI-840A is calibrated a

NOTE: The LI-840A is calibrated at the factory before shipping. It is recommended, however, that you verify the calibration by connecting a gas with known concentration and reading the displayed value. If the instrument needs to be calibrated, follow the steps below.

Connect a dry, CO_2 -free gas to the input air stream. Select **Calibration** from the View menu. Click on the **Zero** tab. Click the **Zero** CO₂. Wait up to 15 minutes for any water vapor to clear from the optical cell, then click the **Zero** H₂O button. After several seconds, a message will appear that indicates the IRGA is zeroed.

Set Conne

Set the analyzer span.

Connect a span gas with known concentration of CO_2 to the input air stream. Choose a span gas that will be below or close to the highest reading expected. Select the **Span** tab. Enter the value of the span gas in ppm. Click on **Span CO₂**. A message will appear indicating when the span calibration is completed.

Connect an airstream with a known water vapor concentration (such as may be generated by the LI-610 Portable Dew Point Generator). Enter the dew point value of the H_2O span gas, in degrees C, and click on **Span H₂O**.

Connect a second span gas with a known concentration of CO_2 to the input air stream. If the reading is not within 1% of the expected value, perform a second span: use a span gas with a concentration that is higher than the highest expected reading. Under the **Span 2** tab enter the 2nd span gas concentration and click **Span CO**₂.

Check the H_2O reading with a water vapor concentration that is higher than that used to set the H_2O span. If the reading is outside the acceptable range, perform a secondary H_2O span. Click **Close**.

12 Set up the logging parameters.

Select **Options** from the Logging menu. Choose the File Header(s) to be output with the data, the values to be logged, the field delimiter, and log frequency. Click **OK**. Choose **Start** from the Logging menu. Enter a file name for the data file. Click **Save**.

13

Set up the charting parameters, if desired.

Select **Charting** from the View menu. Enter the value to be plotted, the Y-Axis min and max values, X-Axis max, and time units to be displayed on each plot. Click **Start** to begin recording data. Graphs with data will appear in the window and data will be logged to the file designated in Step 11 above.

14

Choose Stop from the Logging menu to stop recording when you are done collecting data.

2 Power On

Power On

The LI-840A can be connected to a constant source of AC power using a transformer that provides 12-30VDC (14W maximum) to the LI-840A. If the LI-840A is powered with a voltage below 10.5 volts, the analyzer will not power up, and the Low Battery LED on the top panel will illuminate. The LI-840A will continue to operate with a low battery; there will, however, be a corresponding reduction in performance.

Alarms

The LI-840A is equipped with high and low alarms, which can be configured as open collector or 0-5V output (TTL levels) using jumpers on the main PC board (Figure 2-1 on the next page). The default configuration is 0-5V output. Figure 2-2 on page 2-3 shows the position of the jumpers for each of the two alarm conditions.

The alarms can be used for either CO_2 or H_2O concentrations; only one can be chosen at a time, however.

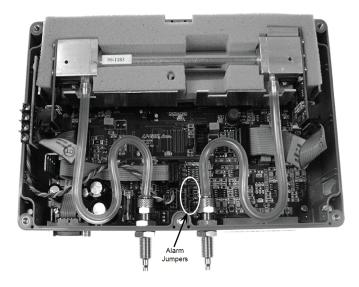


Figure 2-1. Location of alarm jumpers LK1 and LK2.


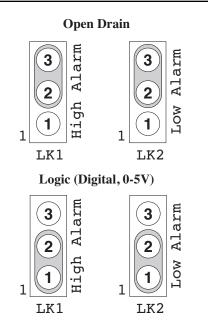


Figure 2-2. Position of jumpers for open drain and 0-5V output.


In addition, a "dead band" value can be set in software for both high and low alarms. To understand how the alarms and dead band values work, look at the diagram below.

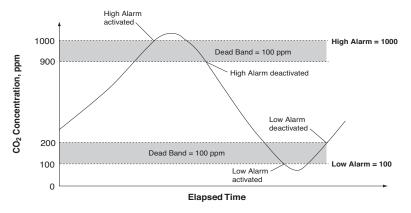

In this chart, the low and high alarm values are set to 100 ppm and 1000 ppm, respectively. The dead band value in both alarms is set to 100 ppm. When the CO_2 concentration reaches 1000 ppm, the high alarm is activated, and remains active until the concentration drops below 900 ppm. When the CO_2 concentration falls below 100 ppm, the low alarm is activated, and remains active until the concentration rises above 200 ppm. Your choice for the dead band value(s) depends on your application and the fluctuation in CO_2 concentrations over time.

Note: Increasing the signal average value in software (see Settings Window; Options; Filter in Section 3) can help reduce fluctuations in readings.

Alarm LEDs can be viewed on the top panel of the LI-840A. Terminals 3 and 5 on the terminal strip are also connected to the High and Low alarms, respectively. This is useful for connecting an audible alarm, for example, or a relay switch to operate another device that will raise or lower the CO_2 concentration to the desired level. The schematic diagram below shows how the high alarm could be connected to a relay switch that triggers an exhaust fan in a greenhouse environment. These relays could also be used to trigger devices such as automatic dialers, alarms, pumps, and valves in industrial and other environments.

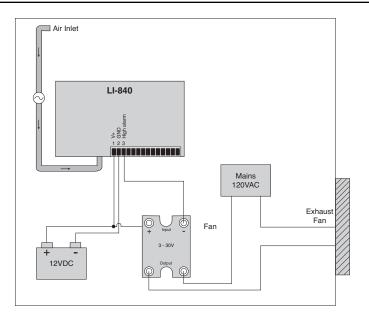

A list of suppliers of electronic relay switches can be found in Appendix D.

NOTE: Consult your local electrical codes before wiring, and/or have a professional electrician wire your application.

3 Operation

Installing the PC Communication Software on Your Computer



The software application is shipped on CD. It requires that your computer have an RS-232 serial (COM) interface, and Windows® XP/Vista/7. The program runs on any PC compatible computer.

An installation menu will start when you insert the CD into your CD-ROM drive. Select *Install Software* from the menu and follow the instructions. If installation does not start, select **Run** from the Windows *Start* menu, and select the LI840Setup.exe file on the CD. When the software has finished the installation procedure, a program icon will be placed in the Programs menu.

NOTE: To remove the software, go to the Control Panel and select *Add/Remove Programs* (XP) or *Uninstall a program* (Vista/7). Choose LI840 from the list of programs and click the Add/Remove button. If your computer has the older LI840 software, be sure to use the correct program when configuring the LI-840A.

Setting the Communication Parameters

LI-840A communication parameters are set automatically when the program is run.

Cabling

The serial cable (null modem) included has 9-pin connectors on both ends; either end plugs into the 9-pin connector (Serial I/O) on the front panel of the LI-840A. Use this cable to interface with your computer's 9-pin serial port. If you want to interface to a computer with a 25-pin serial port, a 9-pin to 25-pin adapter must be used. To connect to a computer with a USB port, a 9-pin serial to USB adapter must be used.

RS-232 Output

Data from the LI-840A can be transferred to a computer for analysis, printing or storage using the RS-232 interface. The LI-840A RS-232 port is configured as Data Terminal Equipment (DTE) with no hardware handshaking, and is bi-directional, meaning information can be transferred both into and out of the LI-840A.

The application software that comes with the LI-840A is used to view data and configure the LI-840A from a PC.

Initial Setup

Click on the LI-840A program icon to start the program. The LI-840A Main Window appears. Select **Connect** from the File menu, or click on the connect icon on the toolbar, beneath the File menu. You are asked to select the serial port to which the LI-840A is connected, and the output interval at which data are output (0.5 to 20 seconds, in 0.5 second increments).

🍯 Connect 📃 🗖 💌
Connection
Com Port: 1
Output Data Every: 0.5 💌 second(s)
Connect Cancel

Choose a COM port and the output interval and click **Connect**. If the instrument is connected properly, data will begin to appear in the Main window (below):

🍯 Ц-840A CO2/H20 Gas Analyzer	
File View Logging Help	
=0= =1 🗭 🗆 🗆 🖬 🖓 🖓	
C02:	H2O:
390.8ppm	14.320ppt
Cell Temperature:	Cell Pressure:
50.2 c	96.3 kPa
Status:	DAC Output
Output Rate: 1 record every 0.5 sec	DAC 1: CO2 (ppm)
Heater: ON	DAC 2: H2O (ppt)
Pressure Comp.: ON	Range: 2.5 v
Alarms: OFF	PC Logging:
Filter: 1 sec	Logging: OFF
	Exit

The Main Window displays the CO₂ concentration (ppm) and H_2O concentration (°C or mmol/mol), cell temperature and pressure, and the status of various LI-840A parameters. There are also three menus used to configure the LI-840A, perform zero and span calibrations, and set up the parameters for recording data.

Note that you can view H_2O concentrations in units of degrees C or mmol/mol. To change the units in the Main window, go to the View menu and toggle the **H2O in ppt** menu item on and off. Alternatively, you can click on the H_2O units being displayed in the Main window to change the display.


Parameter	Description
Cell Temperature	Temperature (°C) in the LI-840A optical cell. This value should remain near 50 °C when the heater is turned ON.
Cell Pressure	Barometric pressure (kPa) measured in the LI-840A optical cell. This value can be used to correct measurements for the effects of atmospheric pressure.
Output Rate	Displays data output rate chosen at connection.
Heater	Shows status of heater (ON/OFF), which is used to maintain the optical cell at a constant 50 °C. The heater is turned ON/OFF in the Settings window.

The LI-840A parameters displayed in the Main window are as follows:

Parameter	Description
Pressure Comp.	Shows whether pressure compensation is ON/OFF. When ON, meas- urements are corrected for the effects of pressure fluctuations on gas density and band broadening.
Alarms	Shows status of High and Low alarms, whose values are set in the Set- tings window. Alarms are enabled (ON), or disabled (OFF).
Filter	Shows the current value for software signal averaging, set in the Settings window. The filter can be set from 0 (no signal averaging) to 20 seconds.
DAC Output	Displays source channels for DAC1 and DAC2, and the voltage output range (0-2.5V or 0-5V).
PC Logging	Shows "Off" if data are not being logged to a file, or "Logging" when data are being logged to a file.

Using the Toolbar

The toolbar in the Main window contains shortcuts for some of the commonly used menu items:

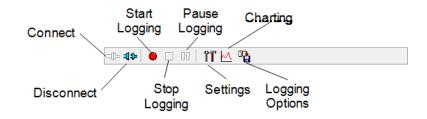

Settings Window - Setting Operational Parameters

The Settings window contains parameters related to initial setup of the LI-840A, including signal filtering, high and low alarm setup, heater and pressure compensation options, and DAC output sources. Choose **Settings** from the View menu (or click on the toolbar icon) to open the Settings window:

Disable the alarms here.	🍯 Settings		Select the DAC output
set high and low alarm values, and high and low	Alarms	DAC Output	 range with these radio buttons.
values, and high and low	Enable Alarms	Range: 0.0v - 2.5v	
Use these radio buttons to choose CO ₂ or H ₂ O as	Source:	DAC 1 Source: H20 ppt	
the alarm source.	High Alarm: 600 ppm	0v:	DAC output sources. Choose data to be
	High Dead Band: 50 ppm	5v:	output to DAC 1 and/or DAC 2.
The heater should be enabled to maintain the	Low Alarm: 400 ppm	DAC 2 Source: CO2 -	
optical bench at a constant 50 °C.	Low Dead Band: 10 ppm	0v: 300	CO ₂ range output by fully scalable DACs.
		5v: 700	Select the desired concentration range
Pressure compensation should be enabled to	Options		that voltage outputs should correspond with.
compensate for gas density changes caused	Enable Heater Fressure Compensation		In this example, 0v
by variations in pressure.	Filter 5 Sec -		represents 300 ppm and 5v is 700 ppm CO ₂
Software signal filtering			The measurement range for CO ₂ with the
is available from 0-20 seconds.		OK Cancel	LI-840A is 0-20,000 ppm.
Seconds.			

Options

Enable Heater

The Heater should be enabled to maintain the optical bench at a constant 50 °C.

Pressure Compensation

Enable the Pressure Compensation check box to automatically correct gas concentration values for changes in cell pressure. In most cases this should be turned on; disabling this feature means that no pressure correction is desired in the gas concentration calculations, which can lead to erroneous measurements.

Filter

Shows the current value for software signal averaging. The filter can be set from 0 (no signal averaging) to 20 seconds. Note that increasing the filter decreases instrument noise but reduces the response time.

DAC Output

Analyzer output for up to 2 values (none, CO_2 , H_2O mmol/mol, H_2O °C, Cell Temperature, or Cell Pressure) can be recorded by connecting a logging device to the terminal strip on the front of the analyzer. Output is linear, and is selectable at 0-5V or 0-2.5V. Choose the full-scale voltage output range and the source for DAC 1 and/or DAC 2. You can also choose a range over which to scale the output(s); you can achieve better resolution by scaling the output(s) over a narrower range when appropriate.

The DACs in the LI-840A are bipolar, and will go slightly negative (\sim -0.100V). This can happen, for example, if the cell becomes contaminated or from small random perturbations when the CO₂ concentration is near zero. See Section 5, Cleaning the Optical Bench for instructions on cleaning the cell should it become contaminated.

Voltage output is measured by attaching the positive lead from the logging device to terminal 9 (V Out 1), or terminal 7 (V Out 2) on the LI-840A terminal strip. Connect the negative lead to position 10 (GND), or position 8 (GND).

The CO_2 or H_2O concentration can be calculated from the DAC output voltage as follows:

$$Concentration = (X_F - X_Z) \frac{V}{V_{\text{max}}} + X_Z$$
3-1

where V is the measured voltage, X_F is the full scale value for CO₂ or H₂O output (entered in the Settings Window as the 2.5 or 5V value, up to 20,000 ppm for CO₂), X_Z is the zero value entered, and Vmax is the full scale DAC output voltage selected (5V or 2.5V).

Example: The DAC output range selected is 0-5V (V_{max}), the zero CO₂ value entered is 0 ppm (X_Z), the full scale CO₂ value entered is 2000 ppm (X_F), and the measured output voltage (V) is 2.9V. To calculate the CO₂ concentration from Equation above,

$$CO_2 = (2000) \frac{2.9}{5} + 0$$
 3-2

= 1160 ppm.

Example: The DAC output range selected is 0-5V (V_{max}), the zero CO₂ value entered is 1000 ppm (X_Z), the full scale CO₂ value entered is 2000 ppm (X_F), and the measured output voltage (V) is 2.9V. To calculate the CO₂ concentration from Equation above,

$$CO_2 = (2000 - 1000) \frac{2.9}{5} - 1000$$
 3-3

= 1580 ppm.

The H₂O concentration (mmol/mol) can be calculated from the DAC output voltage as follows:

$$H_2O = (X_F - X_Z) \frac{V}{V_{\text{max}}} + X_Z$$

$$3-4$$

where V is the measured voltage, X_F is the full scale value for H₂O output (entered in the Settings Window as the 2.5 or 5V value, up to 80 parts per thousand), X_Z is the zero value entered, and V_{max} is the full scale DAC output voltage selected (5V or 2.5V).

Example: The DAC output range selected is 0-5V (V_{max}), the zero H₂O value entered is 0 ppm (X_Z), the full scale H₂O value entered is 80 mmol/mol (X_F), and the measured output voltage (V) is 2.9V. To calculate the H₂O concentration from Equation 3-4 above,

$$H_2O = (80 - 0)\frac{2.9}{5} + 0$$
3-5

= 46.4 mmol/mol.

Example: The DAC output range selected is 0-5V (V_{max}), the zero H₂O value entered is 10 mmol/mol (X_Z), the full scale H₂O value entered is 60 mmol/mol (X_F), and the measured output voltage (V) is 2.9V. To calculate the H₂O concentration from Equation 3-4 above,

$$H_2O = (60 - 10)\frac{2.9}{5} + 10$$
3-6

= 39.0 mmol/mol.

Converting Voltage Output to Cell Temperature

Cell temperature can be calculated from the DAC output voltage as follows:

$$T(^{\circ}C) = (X_F - X_Z)\frac{V}{V_{\text{max}}} + X_Z$$
3-7

where V is the measured voltage, X_F is the full scale value for temperature output (entered in the Settings Window as the 2.5 or 5V value, up to 100 °C), X_Z is the zero value entered, and V_{max} is the full scale DAC output voltage selected (5V or 2.5V).

Converting Voltage Output to Cell Pressure

Cell pressure can be calculated from the DAC output voltage as follows:

$$Pressure(kPa) = (X_F - X_Z) \frac{V}{V_{\text{max}}} + X_Z$$
³⁻⁸

where V is the measured voltage, X_F is the full scale value for cell pressure output (entered in the Settings Window as the 2.5 or 5V value, up to 115 kPa), X_Z is the zero value entered, and Vmax is the full scale DAC output voltage selected (5V or 2.5V).

Converting Current Output to ppm CO₂ or mmol/mol H₂O

Current output can be measured by connecting the positive input of the data logging device to positions 11 or 13 (4-20 mA 2 or 4-20 mA 1), and the negative input to position 12 or 14 (GND). The current output at positions 11 and 13 is non-isolated, and is rated to drive a 250 ohm load.

Note that these channels mirror their respective voltage output channels; current output channel 1 uses the variable chosen for voltage output 1, and current output channel 2 uses the variable chosen for voltage output 2. Similarly, a voltage output of 0 volts corresponds to a current output of 4 mA, and when the voltage outputs are at full scale (2.5 or 5 V), the current output will be 20 mA.

To convert current output (I) to CO_2 or H_2O concentration in your computer or other output device, the following equation can be used:

$$Concentration = (X_F - X_Z) \left(\frac{I-4}{16}\right) + X_Z$$
³⁻⁹

where I is the measured current, X_F is the full scale value for CO_2 or H_2O concentration output (entered in the Settings Window as the 2.5 or 5V full scale value, up to 20,000 ppm for CO_2), and X_Z is the zero value entered.

Example: The zero CO_2 value entered is 0 ppm (X_Z), the full scale CO_2 value entered is 2000 ppm (XF), and the measured current output (I) is 16.25 mA. To convert to ppm CO_2 :

$$CO_2 = (2000) \left(\frac{16.25 - 4}{16}\right) + 0$$
3-10

= 1531.25 ppm.

CO₂ Span Range


The span range entry field determines the maximum CO_2 range over which the LI-840A will make measurements. The factory calibration range for the LI-840A is 0-20,000 ppm. Note that there is a benefit to selecting a narrower range if that range is adequate for your experiment. The narrower range provides better resolution in terms of the digital-to-analog converter (DAC). For example, if you choose a range of 0-1500 ppm, each step of the DAC is able to resolve twice as many data points as compared to the 0-3000 ppm range, which results in better resolution.

Alarms

Allows you to enable/disable the alarms, and enter high, low, and dead band values. A complete discussion can be found in Section 2, Alarms.

Using the Terminal Strip

The terminal strip is located on the front panel of the LI-840A. To connect the wires, insert the bare wire end into the appropriate terminal and tighten the screw above that terminal using the small flat head screwdriver in the spare parts kit. The front face of the terminal strip can be removed to aid in connecting the wires by pulling straight out on the face.

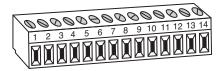


The terminal positions are as follows, reading from left to right:

Terminal	Label	Description	
1	12-30 VDC	Voltage In, 12-30 VDC	
2	GND	Ground	
3	High Alarm	High Alarm	
4	GND	Ground	
5	Low Alarm	Low Alarm	
6	GND	Ground	
7	V OUT 2	Voltage output channel 2	
8	GND	Ground	

Terminal	Label	Description		
9	V OUT 1	Voltage output channel 1		
10	GND	Ground		
11	4-20 mA 2	Current output channel 2		
12	GND	Ground		
13	4-20 mA 1	Current output channel 1		
14	GND	Ground		

Charting Window

Select **Charting** from the View menu to open the Charting window (below). This is the window in which you can set up the parameters for plotting your data. Two charts can be plotted simultaneously, using Y axes on either side of the chart.

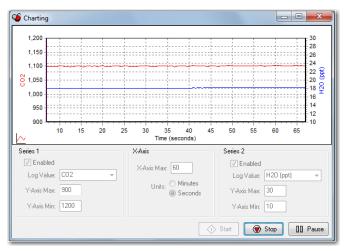

Series 1

The Series 1 options are used to plot a chart with the Y axis on the left side of the chart. Choose the value to be logged and set the maximum and minimum values for the Y axis.

Series 2

The Series 2 options are used to plot a chart with the Y axis on the right side of the chart. Choose the value to be logged and set the maximum and minimum values for the Y axis.


X-Axis Max

Sets the maximum value for the X axis (Time). The units for the X axis can be seconds or minutes.

Press **Start** at any time to view the chart layout and begin displaying data. Note that you must press **Stop** to make changes to the chart parameters, and then press **Start** again to resume data display.

Diagnostic Window

Select **Diagnostics** from the View menu to open the Diagnostics window (below). This window displays the current LI-840A internal software version number, the input voltage, and raw absorption values.

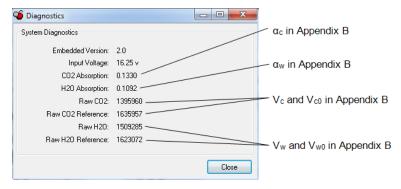

Logging Data

Start Logging

Opens the Log File Destination dialog, where you enter a file name for the data file. The file extension .txt is added automatically.

🍯 Log File Destination				×
- Were the second secon		Search LI840A_2	_0_0c	٩
Organize 🔻 New folder				0
★ Favorites ▲ Name ■ Desktop ▲ ▶ Downloads ▲ ▲ Recent Places ▲	No items match you	Date modified r search.	Туре	
Ubraries ⊡ Documents J Music E Pictures Wideos				
Recomputer	m			
File name: LogFile	m			
Save as type: Text File (*.bxt)				•
Hide Folders		Save	Cancel	

Stop Logging

Stops data logging.


Pause Logging

Pauses logging of data until **Start** is chosen from the Logging menu again, or the Start button on the toolbar is pressed.

Options

Opens the Logging Options window, where you can configure the data output options.

Logging Options	
File Headings	
Date and Time	
V Field Labels	
Log Values	
📝 System Date	
System Time	
CO2	
H20 ppt	
 H20 C Cell Temperature 	
Cell Pressure	
CO2 Absorption	
H20 Absorption	
Field Delimiter	
Space Tab	
 Semicolon 	
Log Frequency	_
Log Rate (seconds): 1	
	OK Cancel

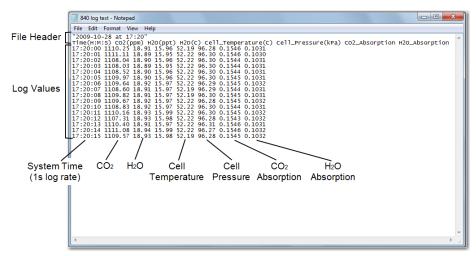

As configured above, the data output would appear similar to that shown below.

Calibration Window - Setting the Zero and Span

Select **Calibration** from the View menu to open the Calibration window. This is the area in which you set the zero and spans of the LI-840A.

Calibration	
Zero Span Span 2 Coefficients Advanced	
Zero CO2	
Last CO2 zero performed on: 11 AUG 2009	
Zero CO2	
Zero H2D	
Last H20 zero performed on: 11 AUG 2009	
Zero H20	
	Close

It is recommended that you perform the zero calibrations first, followed by the span calibrations. To zero, flow a dry, CO_2 -free gas through the LI-840A and make sure the optical cell is completely purged. Press the **Zero CO₂** button under the Zero tab.

The display will show ZERO, and the text in the Calibration window is greyed out. The zero will be set electronically, and the current date will be entered in the "Last CO₂ zero performed on:" field

when completed. Repeat for the H_2O channel. Note: Purge the cell with the H_2O zero gas for up to 15 minutes before zeroing H_2O . Water vapor molecules are "sticky" and it may take up to 15 minutes to reach equilibrium in the optical cell.

Calibration	
Zero Span Span 2 Coefficients Advanced	
Span Last CO2 span performed on:: 11 AUG 2009 Span gas concentration. ppm Span CO2	
Span Last H20 span performed on: 11 AUG 2009	
Span gas concentration: C Span H20	
	Close

To span, connect a span gas of known CO_2 concentration to the input air stream. Make sure the cell is purged. Under the Span tab, enter the value of the span gas and click on **Span CO₂**.

The display will show SPAN, for several seconds and the text in the Calibration window will be greyed out. The span will be set electronically, and the current date will be entered in the "Last CO_2 span performed on:" field when completed. *Note: Ensure that the CO₂ standards are accurate, as this will determine the overall accuracy of the LI-840A.*

Repeat for the H_2O channel using an airstream of known dew point. Note: Purge the cell with the H_2O span gas for up to 15 minutes before spanning H_2O . The LI-610 Portable Dew Point Generator is capable of providing an airstream with a known water vapor concentration.

Two Point Span (Span 2 Tab)

Suppose you have two non-zero calibration gasses (e.g. CO_2 at 300 ppm and 1500 ppm), and you use one of them (say, 1500 ppm) to set the span. If you then flow the other gas through the LI-840A, you would expect it to read within 1% of the expected value (in our example, 300 ±3 ppm). If it does not, you can adjust the calibration for this by performing a Secondary Span, by entering the span gas concentration and clicking the **Span CO₂** (Span 2 tab). After that, the 300 ppm tank should read 300, and the 1500 tank should still read 1500. This can also be performed with H₂O if needed.

Calibration	- • ×
Zero Span Span 2 Coefficients Advanced	
Span	
Last CD2 span 2 performed on: Unknown	
Span gas concentration: ppm	
Span CO2	
Span CO2	
Span	
Last H20 span 2 performed on: Unknown	
Span gas concentration: C	
Span H2D	
Sparrizo	
	Close


In the LI-840A version 2.0 and above, the span value is a linear function of absorptance, and thus has two parameters, a slope and an offset (see Appendix B for details). Normal span setting adjusts the offset and leaves the slope alone. Setting the secondary span adjusts the slope, then adjusts the offset value for the new slope value. The slope and offset values can be viewed on the Advanced page (described below). Span offset values (CO₂ Span and H₂O Span) are typically 1.0 \pm 0.1). Span slope values (CO₂ Span 2 and H₂O Span 2) are typically 0 \pm 0.1). When they are exactly 0, then slope is a constant, and not a function of absorptance.

The CO_2 span slope value is first determined at the factory by setting the span at 200 ppm, and doing an secondary span close to 20000 ppm. This value should remain valid for some time. Things


that are likely to make it change include changing the source and/or changing the detector, a dirty optical cell, and possibly even taking the optical bench apart for cleaning.

Considerations for performing a secondary span:

- Zero first.
- After zeroing, do a (normal) span.
- The span and secondary span concentrations should be as far apart as possible and they should bracket the concentration range of interest to you. (e.g. 200 and 20000, or 300 and 1500, etc.). "Far apart" means in absorptance, not concentration. We recommend absorptance differences of at least 0.1 when choosing concentrations for setting span and secondary span (see the figure below).

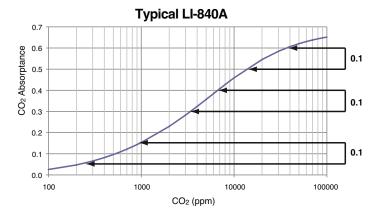

For CO₂, The difference in absorptance between 200 and 1000 ppm is typically about 0.1. There is also a 0.1 absorptance difference between 3500 and 7000 ppm, and between 15,000 ppm and 40,000 ppm. For H₂O, it is a little harder to get a 0.1 absorptance difference. If you want to do a secondary span at room temperature, for example, the two span values would have to be 5 °C or less, and 22 °C, for example.

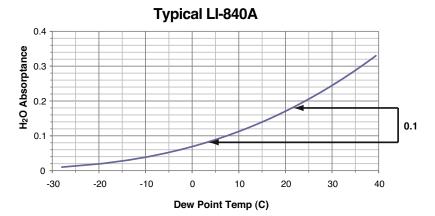

You can do multiple secondary spans after a normal span. For example, span at 200, then do a secondary span at 3000. Then (assuming you didn't like the result for some reason), you can do another secondary span at 2000 ppm. As long as the secondary span is well away (in absorptance) from the previous normal span, it is OK to do as many consecutive secondary spans at various concentrations as you'd like.

Coefficients Page

The LI-840A uses a double rectangular hyperbola for the CO_2 calibration, and a third order polynomial for H_2O calibration. The Coefficients page displays these factory-determined calibration coefficients, as well as those for band broadening and cross sensitivity, and coefficients for the internal pressure transducer. These coefficients are fixed at the factory.

Zero S	òpan	Span 2	Coefficients	Advanced	1				
A3:	5.364 1.040	2600E1 2897E1 1098E2 2500E2			H2O A1: A2: A3:	2.8746101E- 2.8455300 -1.6282000			
Pressu A0: A1:	re				Bano	ctions d Broadening: iss Sensitivity:	1.4500	3	
			Send Coel	ificients		Cancel]		

The calibration coefficients are unique to each instrument, and may be found on the calibration sheet shipped from the factory. In most cases you will not want to alter any of these values.

A description of the band broadening can be found in Appendix B, Equation Summary.

Advanced Page

The calibration constants for CO_2 and H_2O zero and span calibrations are found on the Advanced calibration page. These constants are stored to a file on the computer by clicking on **Save to Defaults**. Click **Load Defaults** to restore the values in this window using the file on the PC. If new constants are entered in this window, click **Send Constants** to send the values to the LI-840A for implementation.

In most cases you should not change the values in this window; see Appendix B, *Equation Summary* for a description of the terms used in this window.

🧯 Calib	oration					- • ×
Zero	Span	Span 2	Coefficients	Advanced		
Cali	ibration Cor	nstants				1
	CO2 Zero:	8.8867	092E-1		Load Defaults	
(CO2 Span:	1.0114	669		Save to Defaults	
CC	02 Span 2:					
	H20 Zero:	1.0032	711			
ŀ	H2O Span:	9.9719	476E-1			
H2	20 Span 2:					
			Send Co	nstants	Cancel	
						Close

Connection the LI-840A to the LI-1400 Datalogger



The following example shows how you can connect the LI-840A to a datalogging device such as the LI-COR Model LI-1400 DataLogger to collect analog data and convert to meaningful CO_2 and/or H₂O values. As mentioned earlier, voltage output from the LI-840A is linear; converting analog data into ppm CO_2 or mmol/mol H₂O simply requires multiplying the raw mV output by a conversion factor, which is dependent upon the selected voltage output range, and the maximum CO_2 or H₂O resolution available (set in software between 0-20,000 ppm). In general, the conversion takes the form:

$$Concentration = (X_F - X_F) \frac{V}{V_{\text{max}}} + X_Z$$
3-11

where V is the measured voltage, X_F is the full scale value for CO₂ or H₂O output (entered in the Settings Window as the 2.5 or 5V value, up to 20,000 ppm for CO₂), X_Z is the zero value entered, and V_{max} is the full scale DAC output voltage selected (5V or 2.5V).

For example, if the voltage output is set for 0-5V (V_{max}), the zero CO₂ value entered is 0 ppm (X_Z) and scaled CO₂ range is 2000 ppm (X_F), and the measured voltage output is 2.9V, the equation would take the form:

$$CO_2 = (2000) \frac{2.9}{5} + 0$$
 3-12

= 1160 ppm.

This general equation can be converted into simple multipliers, based on the two available voltage output ranges and the CO₂ (maximum 20,000 ppm) or H₂O (maximum 100% RH) range. These can be any value, depending on what is chose for the 0V and the full scale 2.5V and 5V DAC output values. Each multiplier can then be entered in the data logging device to convert raw voltage to CO₂ or H₂O values. Table 3-1 lists some examples of values for this multiplier; simply choose your DAC output range in the second row, and then select the LI-840A CO₂ range in the first column (the full scale value minus the zero value); follow across to find the appropriate multiplier. Note that this table assumes that the zero value (X_Z) is zero; if it is not zero, the offset will need to be added onto the final result.

	DAC Output Range				
	0-5V	0-2.5V			
X_F-X_Z	Multiplier (ppm/Volts)				
1000 ppm	200	400			
2000 ppm	400	800			
5000 ppm	1000	2000			
20000 ppm	4000	8000			

Table 3-1. Multipliers for Converting Voltage Output to CO₂ Readings (no offset).

To calculate multipliers:

$$Multiplier = \frac{X_F - X_Z}{V_{\text{max}}}$$
3-13

where X_F is the full scale value for CO_2 or H_2O output, X_Z is the zero or offset value entered, and V_{max} is the full scale DAC output voltage selected (5V or 2.5V).

The LI-1400 can be used to monitor voltage signals up to 2.5 volts with 76 microvolt resolution in real time and convert them into meaningful engineering units shown on the display. Alternatively, the LI-1400 can be configured to automatically log these data over extended periods of time and later transfer the results to a computer for further analysis.

Access to the voltage channels requires the 1400-301 Terminal block. Voltage channels are designated by the letter "V" and sequentially numbered V1-4. The lead from LI-840A terminal #7 or #9 should be attached to one of the LI-1400 terminals labeled V1, V2, V3, or V4, while the lead from terminal #8 or #10 should be attached to one of the LI-1400 terminals labeled \downarrow (ground). See the LI-1400 manual for further details.

The following example shows how you can set up the LI-1400 Data Logger to collect raw voltage output from the LI-840A and convert to CO₂ values.

1. Set the voltage output of the LI-840A for 2.5V output and specify the zero and maximum values for CO₂ as described in Section 3, *Analog Output*.

- **2.** In the LI-1400, configure V1 channel as *General* for CO₂.
- **3.** Enter a description, such as CO₂.
- **4.** Set Math = *Poly(nomial)* and press *Ent(er)*.
- **5.** Set description as desired, a0 = the offset (for scaled measurements), <math>a1 = multiplier from Table 3-1 on page 3-35 (or Equation 3-13), and a2-a5 = 0. When finished, press Esc to return to the main configuration list.
- **6.** Set Oper(ator) = none.
- **7.** Enter a *Label* such as *ppm* for the units.
- **8.** Set Average=1 sec or as desired. CO₂ will now be displayed on channel V1.

To log CO₂ automatically, follow the remaining steps:

- 9. Set Log Routine to the desired log routine.
- **10.** Set Calc=Mean.
- **11.** To capture the minimum and maximum CO_2 values, set *MinMax* accordingly.
- **12.** *TCoef* has no effect when *Calc=Mean*. It is used only when integrating.

4 Theory of Operation

System Overview

The LI-840A CO₂/H₂O Gas Analyzer is an absolute, non-dispersive, infrared (NDIR) gas analyzer based upon a single path, dual wavelength, infrared detection system. The CO₂ and H₂O measurements are a function of the absorption of IR energy as it travels through the optical path. Concentration measurements are based on the difference ratio in the IR absorption between a reference and sample signal. Reference and sample channels measure infrared gas absorption in a single path through the use of narrow band optical filters with appropriately selected bands. The CO₂ sample channel uses an optical filter centered at 4.26 micrometers, corresponding to an absorption band for CO₂. The reference channel for CO₂ has an optical filter centered at 3.95 micrometers, which has no absorption due to CO₂. The H₂O sample channel uses an optical filter centered at 2.595 micrometers, corresponding to an absorption band of H₂O. The reference channel for H₂O has an optical filter centered at 2.35 micrometers, which is a non-absorbing region for H₂O. The instrument uses digital signal processing techniques to determine the temperature and pressure corrected CO_2 concentration based on the optical bench signals using a ratiometric computation. The data are passed through a double rectangular hyperbola (CO_2) or 3rd order polynomial (H_2O) that performs linearization of the detector signal to a mole fraction in air given in µmol CO_2 per mole of air (ppm), and mmol H_2O per mole of air (mmol/mol).

Data output is provided in a digital format through an RS-232 interface that supports connection to an external computer. The instrument comes with a Windows® XP/Vista/7 compatible application for instrument configuration, control, data collection and display. Analog signals are available through a terminal block for collection by a data logger or similar means.

Optical Bench System

The LI-840A CO_2/H_2O Gas Analyzer optical path is a thermostatically controlled IR detection system. The optical bench operation is based upon a broad band IR source and a pyroelectric detector. The source is mounted in a parabolic reflector to collimate the light and increase energy throughput down the optical path to the detector. The reflector and optical path are gold plated to further increase energy transmission. The detectors are pyroelectric devices that operate based on thermal energy received. The narrow band optical filters allow only the wavebands of interest to illuminate the detector, allowing for the determination of CO_2 and H_2O concentrations in the presence of other infrared absorbing gases.

The detectors respond to thermal energy, so it is necessary to precisely regulate the detectors' temperature. This allows for differentiation of thermal gradient noise from the received signals from the optical path. The detection system is shown in Figure 4-1 on the facing page.

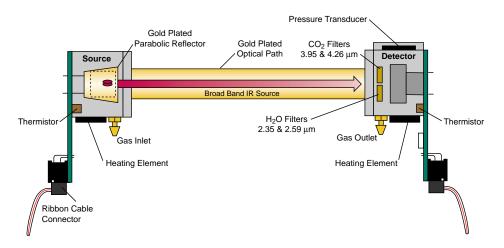


Figure 4-1. Schematic diagram of the LI-840A optical bench.

The optical bench has a thermostat that maintains a constant operating temperature of 50 °C. A feedback loop is used to regulate the optical bench temperature. As shown in Figure 4-1, two thermistors, located in the source and detector housings, measure the present temperature. The thermistors are monitored as part of the control loop to determine corrections necessary in the thermal balance. Two heating elements are the sources of thermal energy into the source and detector housing. The optical path is in mechanical contact with the source and detector housing and thereby achieves thermal equilibrium.

The bench requires approximately 10 minutes to achieve the specified thermal temperature. A longer period of approximately 20 minutes is required to bring the performance of the detection system to within 1 to 2% of reading. As shown in Figure 4-1 on the previous page, the detector housing has a pressure transducer integrated into the housing design. Part of the CO_2 concentration calculation depends on the pressure observed in the optical path, measured with an in-line pressure transducer. Many parameters can affect the pressure and thus the concentration reading. The processing center in the analyzer reads the pressure reading as part of its data collection task and uses this information in the concentration calculation. The gas flow enters the source housing, passes down the optical path and exits at the detector housing. The maximum flow rate for the analyzer is approximately 1 liter/min.

Another key parameter in the concentration calculation is the gas temperature in the optical path. It is assumed in the analyzer operation that the gas temperature will equilibrate to the optical bench temperature (50 °C) by the time it enters the optical path. Since the instrument performs temperature and pressure corrections as part of the concentration calculation, this assumption is very important. To cause the sample air to equilibrate to the optical bench temperature, an airflow pattern is created with sufficient eddy currents to cause thermal equilibration.

The optical bench is mounted in a foam enclosure to accomplish two purposes. The first is to assist in maintaining the controlled thermal environment required for the optical bench as described above. The second function of the foam enclosure is to protect the optical bench from mechanical shock and vibration that might damage the mechanically sensitive components. The foam enclosure is supported by a optical bench "tray" which is in turn attached to the main circuit board (PCB). The optical bench is mechanically de-coupled from the case by creating the s-bend in the Bev-a-line tubing from the external fittings to the optical bench gas inlet and outlet connectors. The electronic interface to the source and detector is achieved through printed circuit boards mounted directly to the source and detector housings, respectively. The PCBs connect the main PCB to the source and detector PCBs via a ribbon cable. This achieves the mechanical isolation desired for the optical bench.


A 14 cm (5.5") optical bench provides a CO_2 measurement range of $0 - 20,000 \mu mol/mol$ (ppm), and an H₂O measurement range of 0-80 mmol/mol.

5 Maintenance


Cleaning the Optical Bench

The LI-840A optical bench can be removed and cleaned if necessary. If the optical path becomes dirty it may become difficult to span the analyzer. Excessive zero drift may also be observed if the optical path becomes dirty. Follow these steps to clean the optical bench:


- **1.** Turn the LI-840A off. Remove the six screws on the LI-840A top panel and remove the cover. Note that these screws are not molded into the case and may fall out.
- **2.** Unscrew the tube retaining nuts on the inner air port fittings. Remove the tubing from both air ports. Leave the tubing connected to the source and detector housings.
- **3.** There are ribbon cables connected to the circuit boards on the source and detector housings, as shown below. Pull straight out on the connector that is attached to each ribbon cable. The optical bench can now be removed from the foam casing.

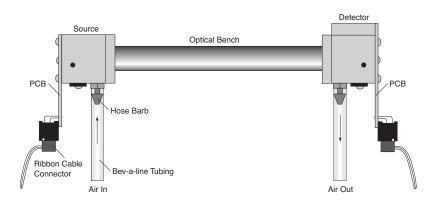

4. There are four screws on the source and detector circuit boards that must be removed. Remove the four screws in the corners of the boards, as shown below (they are slightly larger than the other four screws). Do not remove the remaining four screws.

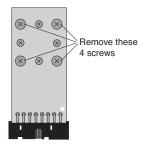


5. The source and detector housings (with attached circuit boards) can now be removed. The bench will appear as shown below. It is a good practice to replace the O-rings when cleaning or replacing the optical path.

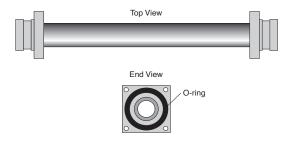

- **6.** There are a number of swabs in the spare parts kit. Dip one end of the swab into a 50:50 ethanol/water solution and carefully swab both ends of the optical bench, until there is no more visible residue. A mild solution of dish washing type soap and water will also work. *Do not use abrasive cleansers, as they can irreparably damage the gold plating on the optical bench.*
- **7.** Use a reflector swab and carefully swab the gold-plated concave surface of the source housing, if necessary.
- **8.** If you need to clean out the hose barbs or replace the tubing connected to the source and detector housings, use a small pair of diagonal cutters to remove the tubing from the hose barbs. Use the cutters to pinch the tubing parallel to the hose barb axis, and then pivot the cutters over the hose barb tip; the tubing will pull off of the hose barb. *Be very careful not to cut the tubing*

or scratch the hose barb with the cutters, as subsequent tubing connections may leak if the hose barb is damaged.


- **9.** Let the optical bench dry. Re-assemble the bench, making sure the O-rings are in place on both ends of the bench. Note that the orientation of the cylinder is not important; either end can be inserted into the source or detector housing.
- **10.** Re-assemble the LI-840A case. Make sure that the foam insulation on the inside top cover is positioned over the optical bench; it is required for thermal stability.
- **11.** Perform zero, span, and secondary span calibrations as described in "*Calibration Window Setting the Zero and Span*" on page 3-25.

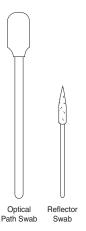

Changing the Fuse

The LI-840A power supply is protected by a 2A 250V, 5×20 mm fast-blow type fuse located inside the case. If the battery fails to power the LI-840A, and will not light the Power LED on the top panel, check to see if the fuse has blown.

To check the fuse, remove the six screws on the top of the LI-840A. The fuse is located on the main circuit board, near the **Flow In** port, as shown below. Replacement fuses (part #439-03952, in the spares kit) plug into the fuse holder; no soldering is required. Replace the fuse and reassemble the LI-840A case.

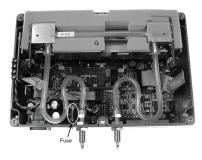

Note on Infrared Source Replacement

The infrared source in the LI-840A is rated for more than 18,000 hours of life. Should the source fail, a replacement can be obtained from LI-COR Biosciences. The source can be replaced in the field; however, it should be noted that changing a source has the potential of degrading the accuracy specification for CO_2 to as much as 3% of reading. The H₂O channel should remain within the published specification.

If the source fails, the instrument can be returned to LI-COR Biosciences for replacement and factory recalibration. This will ensure that the accuracy stays within the factory specifications. Alternatively, if you have two span gasses, you can adjust the calibration yourself. See *"Two Point Span* (Span 2 Tab)" on page 3-27.

Specifications

\mathbf{CO}_2^1

Measurement Range: 0-20,000 ppm Accuracy: Better than 1.5% of reading **Calibration Drift Zero Drift:** <0.15 ppm/°C **Span Drift:** < 0.03 %/°C Total Drift at 370 ppm: <0.4 ppm/°C RMS Noise at 370 ppm with 1 sec signal filtering: <1 ppm Sensitivity to water vapor: < 0.1 ppm CO₂/mmol/mol H₂O H₂O Measurement Range: 0-60 mmol/mol Accuracy: Better than 1.5% of reading **Calibration Drift Drift² at 0 mmol/mol:** <0.003 mmol/mol/°C **Span Drift³ at 10 mmol/mol:** <0.03 %/°C Total Drift⁴ at 10 mmol/mol: <0.009 mmol/mol/°C **RMS Noise at 10 mmol/mol with 1 sec signal filtering:** <0.01 mmol/mol

Sensitivity to CO₂: <0.0001 mmol/mol H₂O/ppm CO₂ Measurement Principle: Non-Dispersive Infrared Traceability: Traceable gases to WMO standards for CO₂. NIST traceable LI-610 Portable Dew Point Generator for H₂O Pressure Compensation Range: 15 kPa-115 kPa Maximum Gas Flow Rate: 1 liter/min **Output Signals:** Two analog voltage (0-2.5V or 0-5V) and two current (4-20mA) Digital: TTL (0-5V) or Open Collector DAC Resolution: 14-bits across user-specified range Source Life: ~18,000 Hours Power Requirements: Input Voltage 12-30 VDC 1.2A @ 12V (14W) maximum during warm-up with heaters on 0.3A @ 12V (3.6W) average after warm-up with heaters on **Operating Temperature Range:** -20°C to +45°C Relative Humidity Range: 0 to 95% RH, Non-Condensing **Dimensions:** $8.75'' \times 6'' \times 3'' (22.23 \times 15.25 \times 7.62 \text{ cm})$ Weight: 2.2 lbs (1 kg)

Internal Optical Cell Volume: 14.5 mL

¹Specifications subject to change without notice
²Zero drift is the change with temperature at 0 concentration.
³Span drift is the residual error after re-zeroing following a temperature change.
⁴Total drift is the change with temperature without re-zeroing or re-spanning.

B Equation Summary

General Form

The LI-840A computes CO₂ and H₂O concentrations using an equation of the form

$$c = f(\alpha g(\alpha, P)S(\alpha)) (T + 273.15)$$
B-1

where c is concentration, f() is the calibration function, α is the absorptance, $g(\alpha, P)$ is the pressure correction, $S(\alpha)$ is the span, and T is the temperature (°C) of the gas in the cell, typically 51.5 °C. Absorptance is computed from

$$\alpha = \left(1 - \frac{V}{V_o}Z\right)$$
B-2

where V and V_o are the raw detector sample and reference readings, and Z is the zeroing parameter. Span is a linear function of absorptance.

$$S(\alpha) = S_0 + S_1 \alpha$$
 B-3

H₂O Equations

Absorptance α_w for water vapor is computed from

$$lpha_w = \left(1 - rac{V_w}{V_{wo}} Z_w
ight)$$
B-4

where V_w and V_{wo} are the sample and reference raw detector readings and Z_w is the zero parameter. The pressure correction for water vapor is an empirical function $g_w()$ of absorptance and pressure P:

$$g_w(\alpha_w, P) = \frac{P_o}{P(1+0.8\alpha_w(\frac{P_o}{P}-1))}$$
B-5

The value of P_o is 99 kPa. When the pressure correction is not enabled, g_w () is simply 1.0. Values for g_w () can be viewed on the diagnosites screen (Section 3).

Water vapor concentration W (mmol mol⁻¹) is computed from

$$W = f_w \left(lpha_w g_w(lpha_w, P) S_W \left(lpha_w
ight)
ight) (T + 273 - 15)$$
 B-6

where $f_w(x)$ is a third order polynomial whose coefficients are given on the calibration sheet.

$$f_w\left(x
ight) = a_{w1}x + a_{w2}x^2 + a_{w3}x^3$$
 B-7

CO₂ Equations

The measurement of CO_2 is a bit more complicated than for H_2O because of the influence of water vapor. There is a slight direct cross sensitivity in the CO_2 signal to H_2O . This is measured at the factory and accounted for in the computation of absorptance (Equation B-8). There is also a band broadening effect that is accounted for in the computation of concentration (Equation B-12).

 CO_2 absorptance α_c is computed from

$$\alpha_{c} = \left(1 - \left(\frac{V_{c}}{V_{co}} + X_{wc}\left(1 - \frac{V_{w}}{V_{wo}}\right)\right)Z_{c}\right)$$
B-8

where V_c and V_{co} are the raw detector signals for sample and reference, Z_c is the CO₂ zero parameter, and X_{wc} is a cross sensitivity parameter for the effect of water vapor on CO₂. Its value is reported on the calibration sheet as "XS=".

The empirical pressure correction function g_c () depends on CO₂ absorptance and pressure:

When $P = P_o, g_c() = 1$.

When $P < P_o$

$$g_c\left(\alpha_c,P\right)=X$$

$$X = \frac{1}{A + B\left(\frac{1}{z - \alpha_c} - \frac{1}{z}\right)} + 1$$

$$A = \frac{1}{a(p-1)}$$
B-9

$$B=rac{1}{rac{1}{b+cp}+d}$$
 $p=rac{P_0}{P}$

where a = 1.10158, $b = -6.1217 * 10^{-3}$, c = -0.266278, d = 3.69895, and z is the asymptotic value of absorptance, obtained from the calibration coefficients (see Equation B-13).

$$z = a_{c1} + a_{c3}$$
B-10

When
$$P > P_o$$

 $g_c(\alpha_c, P) = \frac{1}{X}$
 $P = \frac{P}{P_0}$
B-11

where X, A, and B are computed as in Equation B-9. The variable g_c () is viewable on the Diagnostics screen (see " Operation" on page 3-1).

 CO_2 concentration C (µmol mol⁻¹) is computed from

$$C = f_c \left(\frac{\alpha_c g_c(\alpha_c, P)}{\psi(W)} S_c(\alpha_c) \right) \psi(W) (T + 273.15)$$
B-12

where $f_c(x)$ is a function whose inverse is a double rectangular hyperbola, and whose coefficients (a1...a4) are given on the calibration sheet.

$$f_c^{-1}(C) = \frac{a_{c1}C}{a_{c2}+C} + \frac{a_{c3}C}{a_{c4}+C}$$
B-13

Solving Equation B-13 for C yields the calibration function

$$f_c\left(x\right) = \frac{\left(a_2a_3 + a_1a_4\right) - \left(a_2 + a_4\right)x - \sqrt{\left(a_2 - a_4\right)^2 x^2 + Dx + \left(a_2a_3 + a_1a_4\right)^2}}{2(x - a_1 - a_3)}$$
B-14

Where

$$D = 2(a_2 - a_4)(a_1a_4 - a_2a_3)$$
B-15


 $\psi(W)$ accounts for band broadening by water vapor, and is viewable on the Diagnostic screen (Section 3).

$$\psi(W) = 1 + (h(\alpha_c) - 1) \frac{W}{1000}$$
 B-16

The band broadening coefficient $h(\alpha_c)$ has been determined to be 1.45 for the LI-840A for CO₂ concentrations near ambient. At higher concentrations, the value decreases. We capture this behavior with an empirical relationship (Equation B-17).

$$h(\alpha_c) = \frac{1}{(0.64b_w - 0.64)e^{-3\left(\frac{z}{\alpha_c} - 1\right)} + \frac{1}{b_w}}$$
B-17

Where z is from Equation B-10, and b_w is the low concentration band broadening coefficient: 1.45. This is the value shown on the calibration sheet as BB=1.45. The present value of $h(\alpha_c)$ is viewable on the Diagnostics screen (Section 3). The typical relationship between $h(\alpha_c)$ and CO₂ concentration is illustrated below. ('Typical' because the exact relationship depends on the relationship between absorptance and CO₂, which is the calibration curve.)

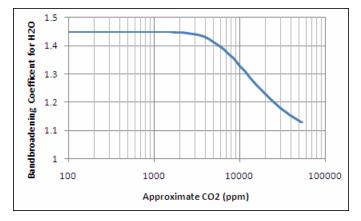

<u>Implementation Note:</u> We formulated Equation B-17 with $0.64b_w - 0.64$ instead of the simple equivalent (0.29) because this allows band broadening corrections to be turned off by setting b_w to 1.

When $b_w = 1$, $h(\alpha_c) = 1$ everywhere. Also, to avoid computational problems (underflows, overflows, and division by zero) we constrain the argument α_c when computing $h(\alpha_c)$ to be

$$0.1 < \alpha_c \le z$$
 B-18

 $\alpha_c - 0.1$ is typically equivalent to about 600 ppm.

Zeroing H₂O

When the command for zeroing water is received, the LI-840A computes the water zero from Equation B-19, where \overline{V}_{wo} and \overline{V}_{wo} are averaged for 5 seconds.

$$Z_w = \frac{\overline{V}_{wo}}{\overline{V}_w}$$
B-19

Zeroing CO₂

When the command for zeroing CO₂ is received, the LI-840A computes the CO₂ zero term from Equation B-20, where \overline{V}_c , \overline{V}_{co} , \overline{V}_{w} , and \overline{V}_{wo} are averaged for 5 seconds.

$$Z_{c} = \frac{1}{\left(\frac{\overline{V}_{c}}{\overline{V}_{co}} + X_{wc} \left(1 - \frac{\overline{V}_{w}}{\overline{V}_{wo}} Z_{w}\right)\right)}$$
B-20

Spanning H₂O

When the command for setting the span for H₂O is received, along with the target concentration W_{T_2} the LI-840A computes S_{w0} from Equation B-21, where $\overline{\alpha}_w$ is averaged over five seconds.

$$S_{w0} = rac{eta_w}{\overline{lpha}_w} - S_{w1}\overline{lpha}_w$$
 B-21

where

1

$$\beta_w = \frac{f_w^{-1}\left(\frac{W_T}{T+273.15}\right)}{g_w(\overline{\alpha}_w, P)}$$
B-22

The instrument retains the following values, which are used for subsequent secondary spans:

$$\alpha_{w1} = \overline{\alpha}_w$$
B-23

$$\beta_{w1} = \beta_w$$
 B-24

Secondary Span H₂O

When the secondary span command for H₂O is received, the instrument computes new values for both S_{w0} and S_{w1} . First, it measures a new $\overline{\alpha}_{w}$ and computes a new β_{w} from Equation B-22. Then, it uses these plus the retained values (α_{w1} and β_{w1} from the previous normal span) to compute

$$S_{w1} = \frac{\frac{\beta_w}{\overline{\alpha}_w} - \frac{\beta_{w1}}{\overline{\alpha}_{w1}}}{\overline{\alpha}_w - \alpha_{w1}}$$
B-25

Given the new span slope S_{wl} , update the span offset S_{w0} by Equation B-21.

Spanning CO₂

When the command for setting the span for CO₂ is received, along with the target concentration C_T , the LI-840A computes S_{c0} from Equation B-26, where $\overline{\alpha}_c$ and \overline{W} are averaged for 5 seconds.

$$S_{c0} = \frac{\beta_c}{\overline{\alpha}_c} - S_{c1}\overline{\alpha}_c$$
 B-26

where

$$\beta_{c} = \frac{f_{c}^{-1} \left(\frac{C_{T}}{(T+273.15)\psi(\overline{W})} \right) \psi(\overline{W})}{g_{c}(\overline{\alpha}_{c}, P)}$$
B-27

Note that

$$\psi(\overline{W}) = 1 + (h(\overline{\alpha}_c) - 1) \frac{\overline{W}}{1000}$$
$$= \left(1 + \left(\frac{1}{(0.64b_w + 0.64)e^{-3\left(\frac{z}{\overline{\alpha}_c} - 1\right)} + \frac{1}{b_w}}\right) \frac{\overline{W}}{1000}\right)$$
B-28

The instrument retains the following values which are used for subsequent secondary spans, if necessary:

$$\alpha_{c1} = \overline{\alpha}_c$$
 B-29

$$\beta_{c1} = \beta_c$$
 B-30

Secondary Span CO₂

When the secondary span command for CO₂ is received, the instrument computes new values for both S_{c0} and S_{c1} . First, it measures a new $\overline{\alpha}_c$ and computes a new β_c from Equation B-27. Then it uses these, plus the retained values (α_{c1} and β_{c1} from the previous normal span) to compute

$$S_{c1} = \frac{\frac{\beta_c}{\overline{\alpha}_c} - \frac{\beta_{c1}}{\alpha_{c1}}}{\overline{\alpha}_c - \alpha_{c1}}$$
B-31

Given the new span slope S_{cl} , update the span offset S_{c0} by Equation B-26.

Symbol Summary

Many of the quantities described in the above section are available from the LI-840A. These are summarized in Table B-1 below.

 Table B-1. Symbol summary


Symbol	XML Tag: <li840>+</li840>	Description
$\alpha_{\rm W}$	<data><h2oabs></h2oabs></data>	H ₂ O absorptance
$g_W(\alpha_w,P)$	<auxdata><pha></pha></auxdata>	Pressure corrected absorptance of H ₂ O
W	<data><h2o></h2o></data>	$H_2O \pmod{mol^{-1}}$
Р	<data><cellpress></cellpress></data>	Pressure (kPa)
Т	<data><celltemp></celltemp></data>	Temp (C)

 Table B-1. Symbol summary (...continued)

Symbol	XML Tag: <li840>+</li840>	Description
α_c	<data><co2abs< td=""><td>CO₂ absorptance</td></co2abs<></data>	CO ₂ absorptance
$g_c(\alpha_c, P)$	<auxdata><pca></pca></auxdata>	Pressure corrected absorptance of CO ₂
С	<data><co2></co2></data>	$CO_2 \ (\mu mol \ mol^{-1})$
V_{w}	<data><raw><h2o></h2o></raw></data>	Raw detector readings
V_{wo}	<data><raw><h2oref></h2oref></raw></data>	
Vc	<data><raw><co2></co2></raw></data>	
V_{co}	<data><raw><co2ref></co2ref></raw></data>	
$a_{w1}a_{w3}$	<poly><h2o></h2o></poly>	Calibration coefficients
$a_{c1}a_{c4}$	<poly><co2></co2></poly>	Calibration coefficients
$Z_{\rm w}$	<cal><h2okzero></h2okzero></cal>	
S_{w0}	<cal><h2okspan></h2okspan></cal>	
\mathbf{S}_{w1}	<cal><h2okspan2></h2okspan2></cal>	
Zc	<cal><co2kzero></co2kzero></cal>	Zero and Span values
Sc0	<cal><co2kspan></co2kspan></cal>	
S _{c1}	<cal><co2kspan2></co2kspan2></cal>	
X _{wc}	<poly><xs></xs></poly>	Cross sensitivity coefficient
b_{w}	<poly><bb></bb></poly>	Band broadening coefficient for low CO ₂
$h(\alpha_c)$	<auxdata><bb_eff></bb_eff></auxdata>	Computed band broadening coeff for the current absorptance
ψ(W)	<auxdata><psi></psi></auxdata>	Band broadening correction factor for water vapor

C Pin Assignments

DB-9 Connector

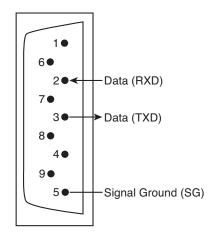

Suppliers

Although the LI-840A has no internal chemicals to replace, the following suppliers may be useful for obtaining chemicals for generating a chemical zero gas. The company names, addresses, and phone numbers are the most current we have at the time of this printing. In some cases the information may change without notice.

Soda Lime (6-12 mesh) and Magnesium Perchlorate (Anhydrous) Mg(ClO₄)₂

GFS Chemicals

www.gfschemicals.com	Soda Lime: Part #66352
P.O. Box 245	Mg(ClO ₄) ₂ : Part #49001 (500g, <8% water)
Powell, OH 43065	
Phone: 614-881-5501	
FAX: 614-881-5989	
Toll free: 800-858-9682	

Fisher Scientific

www.fishersci.com 711 Forbes Avenue Pittsburgh, PA 15219-4785 Phone: 201-467-6400 FAX: 201-379-7415 Toll free: 800-776-7000 Toll free FAX: 800-926-1166

Soda Lime: Part #S201-212 (LI-COR Part #9960-071) **Mg(ClO₄)₂:** Part #M54-500 (500g)

Thomas Scientific

www.thomassci.com P.O. Box 99 Swedesboro, NJ 08085-6099 Phone: 609-467-2000 FAX: 609-467-3087 Toll free: 800-345-2100 Toll free FAX: 800-345-5232

Soda Lime: Part #C703-B76 **Mg(ClO₄)₂:** Part #C260-M61 (Dehydrite, 500g)

Ascarite II

GFS Chemicals	Thomas Scientific	
www.gfschemicals.com	www.thomassci.com	
P.O. Box 245	P.O. Box 99	
Powell, OH 3065	Swedesboro, NJ 08085-6099	
Phone: 614-881-5501	Phone: 609-467-2000	
FAX: 614-881-5989	FAX: 609-467-3087	
Toll free: 800-858-9682	Toll free: 800-345-2100	
	Toll free FAX: 800-345-5232	

P.W. Perkins Co., Inc.	Fisher Scientific	
www.pwperkins.com	www.fishersci.com	
856-769-3525	711Forbes Avenue	
	Pittsburgh, PA 15219-4785	
	Phone: 201-467-6400	
	FAX: 201-379-7415	
	Toll free: 800-776-7000	
	Toll free FAX: 800-926-1166	

Electronic Relay Switches

Crydom Inc.	Potter & Brumfield Products Div.
9525 Chesapeake Dr.	Siemens Electromechanical Components, Inc.
San Diego, CA 92123	200 S. Richland Creek Dr.
800-827-9366	Princeton, IN 47671-0001
FAX: 619-715-7280	info@ae.sec.siemens.com

E

Configuration Grammar

Introduction

The LI-840A communicates exclusively through a serial interface. The following discussion describes how to implement a synchronized communication protocol.

LI-840A Communications and XML

The configuration grammar used to communicate with the LI-840A is based upon a subset of the eXtensible Markup Language (XML). XML relies on the use of tags to "Markup", or give structural rules to a set of data.

A tag is a descriptive identifier, enclosed between a less than (\leq) and a greater than (\geq) symbol, used in part to describe a piece of data. For example, <NAME> is a tag that describes a person's name. Each tag must have a corresponding end tag, denoted by '/'. Extending the example above, the end tag of <NAME> is </NAME>.

Elements are the basic unit of XML content. An element consists of a start tag and an end tag, and everything in between. For example, consider the following element:

<NAME>George</NAME>.

In this example, <NAME> (start tag) and </NAME> (end tag) comprise the markup, and "George" comprises the data. Because XML is extensible, tags can be defined specifically for the data they are meant to describe.

Elements can also contain other elements other than data.

```
<NAME>
<FIRST>George</FIRST>
<LAST>Smith</LAST>
</NAME>
```

In this example, the outermost element <NAME> encompasses two other elements that contain data. All elements combined make up the XML document.

Connecting and Configuring Data Output

The LI-840A communicates through a serial port on the front of the instrument. This port is configured as follows:

Baud Rate: 9600 bps Data Bits: 8 Parity: None Stop Bits: 1

Flow Control: None

After a serial connection is established, the LI-840A will immediately send data out the serial port in the manner in which it was configured previously. In order to reconfigure the LI-840A to output specific data values, you must send the RS-232 portion of the XML grammar to the instrument with the desired values "turned on". To "turn on" the data value, set the value of the element to TRUE. As an example, the following string is sent to the instrument after a connection has been made between the computer (using the Windows® application software) and the LI-840A:

```
<LI840>
<CFG>
<OUTRATE>1</OUTRATE>
</CFG>
<RS232>
<STRIP>FALSE</STRIP>
<ECHO>FALSE</ECHO>
<CELL TEMP>TRUE</CELL TEMP>
<CO2>TRUE</CO2>
```

```
<CO2ABS>FALSE</CO2ABS>
<CELLPRES>TRUE</CELLPRES>
<IVOLT>TRUE</IVOLT>
</RS232>
</LI840>
```

Sending Data to the LI-840A

To send data to the LI-840A, each string must end with a '\n' to ensure that the LI-840A can parse consecutive commands.

After data have been sent to the LI-840A, the instrument replies with:

<LI840><ACK>TRUE</ACK></LI840>

if the XML was received and parsed correctly. If there was an error in the XML, the following is sent:

<LI840><ACK>FALSE</ACK></LI840>.

For example, suppose that you would like to reconfigure the LI-840A to stop outputting cell temperature. Here is the command to send to the LI-840A:

<LI840><RS232><CELLTEMP>FALSE</CELLTEMP></RS232></LI840>

If the command was received correctly, the LI-840A replies with

E-4 Appendix E. Configuration Grammar

<LI840><ACK>TRUE</ACK></LI840>.

Reading Data From the LI-840A

The LI-840A can send data continuously. To determine where one message ends and the next begins, each XML document sent from the LI-840A is delimited with a n' (0x10) character.

Polling the Current State of the LI-840A

The LI-840A can be polled for individual sets of data by sending an XML document with a '?' in place of the set of elements requested. The element sets that can be requested include the data set, the current configuration, and the entire state of the instrument.

Sending this command:

```
<LI840><DATA>?</DATA></LI840>
```

instructs the LI-840A to send the most recent set of data values (as configured) as an XML document.

Sending this command:

```
<LI840><CFG>?</CFG></LI840>
```

instructs the LI-840A to send an XML document containing all of the configuration information, including heater status, filter settings, DACs, and alarms.

To receive the entire state of the instrument as an XML document, send this command: <LI840>?</LI840>.

Calibration

The LI-840A CO_2 calibration (zero and span) can be performed using XML grammar. This is accomplished in three steps:

1. Send the calibration command to the LI-840A.

```
<LI840>
<CAL>
<DATE>{iso date}</DATE>
<CO2SPAN>{int}</CO2SPAN>
<CO2SPAN2>{int}</CO2SPAN2>
<CO2KZERO>{bool}</CO2KZERO>
<co2kspan>{float}</co2Kspan>
<co2kspan2>{float}</co2Kspan2>
</CAL>
</LI840>
```

2. An acknowledgement is returned from the LI-840A if the command was accepted.

```
<LI840><ACK>TRUE</ACK></LI840>
```

3. After a few seconds, all of the calibration information is returned from the LI-840A, indicating that it has finished the zero and/or span.

```
<LI840>
<CAL>
<CO2LASTZERO>{iso date}</CO2LASTZERO>
<CO2KZERO>{float}</CO2KZERO>
<CO2LASTSPAN>{iso date}</CO2LASTSPAN>
<CO2LASTSPAN2>{iso date}</CO2LASTSPAN2>
<CO2KSPAN2{float}</CO2KSPAN>
<CO2KSPAN2>{float}</CO2KSPAN2>
</CAL>
</LI840>
```

If the calibration cannot be performed, an ERROR is sent:

```
<LI840><ERROR>{Error Text}</ERROR></LI840>.
```

To Zero the LI-840A

1. Send the XML command to initiate the zero.

```
<LI840>
<CAL>
<DATE>YYYY-MM-DD</DATE>
<CO2ZERO>TRUE</CO2ZERO>
</CAL>
</LI840>
```

- **2.** Wait for the acknowledgement.
- **3.** Wait for the date to be returned to verify the zero operation succeeded. If the operation fails an <ERROR> will be sent.

To Span the LI-840A

1. Send the XML command to initiate the span.

```
<LI840>
<CAL>
<DATE>YYYY-MM-DD</DATE>
<CO2SPAN>Gas Concentration</CO2SPAN>
</CAL>
</LI840>
```

- 2. Wait for the acknowledgement.
- **3.** Wait for the date to be returned to verify the span operation succeeded. If the operation fails an <ERROR> will be sent.

To Perform a Two Point Span

1. Send the XML command to initiate the first span.

```
<LI840>
<CAL>
<DATE>YYYY-MM-DD</DATE>
```

```
<CO2SPAN>Gas Concentration</CO2SPAN>
</CAL>
</LI840>
```

- **2.** Wait for the acknowledgement.
- **3.** Wait for the date to be returned to verify the span operation succeeded. If the operation fails an <ERROR> will be sent.
- **4.** Send the XML command to initiate the second span.

```
<LI840>
<CAL>
<DATE>YYYY-MM-DD</DATE>
<CO2SPAN2>Gas Concentration</CO2SPAN2>
</CAL>
</LI840>
```

- **5.** Wait for the acknowledgment.
- **6.** Wait for the date to be returned to verify the span operation succeeded. If the operation fails an <ERROR> will be sent.

LI-840A XML Grammar and Element Description

Data Types in the XML Grammar

{val | val |...} The value will be a member of the specified set. | = or. {bool}

Boolean values, TRUE | FALSE.

 $\{float\}$

Floating point values in decimal or exponential notation.

{int}

Integers.

{iso date}

A date in the ISO format. 4 digit year - 2 digit month - 2 digit day. Example: 2002-04-27.

XML Grammar

```
<LI840>
<ACK>{bool}</ACK>
<VER>{string}</VER>
<DATA>
<CELLTEMP>{float}</CELLTEMP>
<CELLPRES>{float}</CELLPRES>
<CO2>{float}</CO2>
<CO2ABS>{float}</CO2ABS>
<H2O>{float}</H2O>
<H2OABS>{float}</H2OABS>
```

Appendix E

```
<H2ODEWPOINT> { float } </H2ODEWPOINT>
     <IVOLT>{float}</IVOLT>
      <RAW>
        <CO2>{int}</CO2>
        <CO2REF>{int}</CO2REF>
        <H2O>{int}</H2O>
        <H2OREF>{int}</H2OREF>
     </RAW>
  </DATA>
<AUXDATA>
  <PCA>{float}</PCA>
  <PHA>{float}</PHA>
  <BB EFF>{float}</BB EFF>
  <PSI>{float}</PSI>
</AUXDATA>
<RS232>
  <CO2>{bool}</CO2>
  <H2O>{float}</H2O>
  <CELLTEMP> {bool } </CELLTEMP>
  <CELLPRES>{bool}</CELLPRES>
  <IVOLT>{bool}</IVOLT>
  <CO2ABS>{bool}</CO2ABS>
  <H2OABS>{bool}</H2OABS>
  <H2ODEWPOINT> {bool } </H2ODEWPOINT>
  <RAW>
     <CO2>{bool}</CO2>
     <CO2REF>{bool}</CO2REF>
     <H2O>{bool}</H2O>
     <H2OREF>{bool}</H2OREF>
```

</RAW> <ECHO>{bool}</ECHO> <STRIP>{bool}</STRIP> </RS232> <CFG> <HEATER> {bool } </HEATER> <PCOMP>{bool}</PCOMP> <FILTER>{int}</FILTER> <OUTRATE>{float}</OUTRATE> <ALARMS> <SOURCE>H20 | CO2<SOURCE> <ENABLED> {bool } </ENABLED> <HIGH>{float}</HIGH> <HDEAD>{float}</HDEAD> <LOW>{float}</LOW> <LDEAD>{float}</LDEAD> </ALARMS> <BENCH> { 5 | 14 } </BENCH> {int} <DACS> <RANGE>{2.5|5.0}</RANGE> <SET1>{float}</SET1> <SET2>{float}</SET2> <D1>{NONE | CO2 | H20 | H20DP | CELLTEMP | CELLPRES }</D1> <D1 0>{float}</D1 0> <D1 F>{float}</D1 F> <D2>{NONE | CO2 | H20 | H20DP | CELLTEMP | CELLPRES } </D2> <D2 0>{float}</D2 0>

```
<D2_0>{float}</D2_F>
```

```
</DACS>
</CFG>
<CAL>
  <DATE>{serial number and iso date}</DATE>
  <CO2ZERO>{bool}</CO2ZERO>
  <CO2SPAN>{int}</CO2SPAN>
  <co2span2>{int}</co2span2>
  <H2OZERO>{bool}</H2OZERO>
  <H2OSPAN>{int}</H2OSPAN>
  <H2OSPAN2>{int}</H2OSPAN2>
  <CO2KZERO>{float}</CO2KZERO>
  <CO2KSPAN>{float}</CO2KSPAN>
  <co2kspan2>{float}</co2span2>
  <H2OZERO>{float}</H2OZERO>
  <H2OKSPAN>{float}</H2OKSPAN>
  <h2okspan2>{float}</h2ospan2>
```

or

```
<CO2LASTZERO>{iso date}</CO2LASTZERO>
<CO2LASTSPAN>{iso date}</CO2LASTSPAN>
<H2OLASTZERO>{iso date}</H2OLASTZERO>
<H2OLASTSPAN>{iso date}</H2OLASTSPAN>
</CAL>
<POLY>
<DATE>{string}</DATE>
<BB>{float}</BB>
<XS>{float}</XS>
<RESET>{bool}</RESET>
```

```
<C02>
        <A1>{float}</A1>
        <A2>{float}</A2>
        <A3>{float}</A3>
        <A4>{float}</A4>
     </C02>
     <H2O>
        <A1>{float}</A1>
        <A2>{float}</A2>
        <A3>{float}</A3>
     </H2O>
     <PRESS>
        <A0>{float}</A0>
        <A1>{float}</A1>
     </PRESS>
  </POLY>
   <ERROR>{string}</ERROR>
</LI840>
```

Tag	Parent	Value(s)	R/W	Comments
<li840></li840>	N/A (root)	<ack></ack>		<li840> is the root tag for all XML statements.</li840>
		<ver></ver>		A ? outputs the entire XML document.
		<error></error>		
		<data></data>		
		<auxdata></auxdata>		
		<rs232></rs232>		
		<cfg></cfg>		
		<cal></cal>		
		<poly></poly>		
		?		

<LI840>?</LI840> <LI840><ACK>TRUE</ACK></LI840>

Тад	Parent	Value(s)	R/W	Comments
<data></data>	<li840></li840>	<co2></co2>		<data> contains all data values sent from the</data>
		<co2abs></co2abs>		LI-840A. All of the elements within the
		<h2o></h2o>		DATA tag are readable only. A ? requests all
		<h2odewpoint></h2odewpoint>		data values to be output.
		<h2oabs></h2oabs>		
		<celltemp></celltemp>		
		<cellpres></cellpres>		
		<ivolt></ivolt>		
		<raw></raw>		
		?		

<L1840><DATA><CO2>2.34e2</CO2><IVOLT>1.5e2</IVOLT></DATA></L1840><L1840><DATA>?</DATA></L1840>

Тад	Parent	Value(s)	R/W	Comments
<co2></co2>	<data></data>	Float	R	CO ₂ in ppm
<co2abs></co2abs>	<data></data>	Float	R	CO ₂ absorption
<h<sub>2O></h<sub>	<data></data>	Float	R	H ₂ O in mmol/mol
<h2odewpoint></h2odewpoint>	<data></data>	Float	R	H ₂ O in °C
<h2oabs></h2oabs>	<data></data>	Float	R	H ₂ O absorption
<celltemp></celltemp>	<data></data>	Float	R	Cell temperature
<cellpres></cellpres>	<data></data>	Float	R	Cell pressure
<ivolt></ivolt>	<data></data>	Float	R	Input voltage
<raw></raw>	<data></data>	<co2></co2>		Raw detector readings
		<co2ref></co2ref>		
		<h2o></h2o>		
		<h2oref></h2oref>		

<LI840><DATA><CELLTEMP>5.16E1</CELLTEMP><CELLPRES>9.742E1</CELLPRES><CO2>6.17E2</CO2><CO2ABS>8.94E2</CO2ABS>DATA></LI840>

Тад	Parent	Value(s)	R/W	Comments
<co2></co2>	<raw></raw>	Integer	R	
<co2ref></co2ref>	<raw></raw>	Integer	R	
<h2o></h2o>	<raw></raw>	Integer	R	
<h2oref></h2oref>	<raw></raw>	Integer	R	

Тад	Parent	Value(s)	R/W	Comments
<auxdata></auxdata>	<li840></li840>	?		<auxdata> contains all auxiliary data values that can be retrieved from the LI-840A. All of the elements witin the AUXDATA tag are readable only. A ? requests all data values to be output.</auxdata>

Тад	Parent	Value(s)	R/W	Comments
<pca></pca>	<auxdata></auxdata>	Float	R	
<pha></pha>	<auxdata></auxdata>	Float	R	
<psi></psi>	<auxdata></auxdata>	Float	R	
<bb_eff></bb_eff>	<auxdata></auxdata>	Float	R	

Тад	Parent	Value(s)	R/W	Comments
<rs232></rs232>	<li840></li840>	<co2></co2>		Setting <rs232> values will determine</rs232>
		<co2abs></co2abs>		what values are output in <data></data>
		<h2o><h2odewpoint></h2odewpoint></h2o>		
		<h2oabs></h2oabs>		
		<celltemp></celltemp>		
		<cellpres></cellpres>		
		<ivolt></ivolt>		
		<raw></raw>		
		<echo></echo>		
		<strip></strip>		

Тад	Parent	Value(s)	R/W	Comments
<co2></co2>	<rs232></rs232>	TRUE FALSE	R/W	CO ₂ in ppm
<co2abs></co2abs>	<rs232></rs232>	TRUE FALSE	R/W	CO ₂ absorption
<h2o></h2o>	<rs232></rs232>	TRUE FALSE	R/W	H ₂ O in mmol/mol
<h2odewpoint></h2odewpoint>	<rs232></rs232>	TRUE FALSE	R/W	H ₂ O dew point
<h2oabs></h2oabs>	<rs232></rs232>	TRUE FALSE	R/W	H ₂ O absorption
<celltemp></celltemp>	<rs232></rs232>	TRUE FALSE	R/W	Cell temperature
<cellpres></cellpres>	<rs232></rs232>	TRUE FALSE	R/W	Cell pressure
<ivolt></ivolt>	<rs232></rs232>	TRUE FALSE	R/W	Input voltage
<raw></raw>	<rs232></rs232>	<co2></co2>		Raw detector readings
		<co2ref></co2ref>		
		<h2o></h2o>		
		<h2oref></h2oref>		
<echo></echo>	<rs232></rs232>	TRUE FALSE	R/W	Echo commands sent to LI-840
<strip></strip>	<rs232></rs232>	TRUE FALSE	R/W	Strip XML from all data sent

<LI840><CFG><OUTRATE>0.5</OUTRATE></CFG><RS232><STRIP>FALSE</STRIP> <ECHO>FALSE</ECHO><CELLTEMP>TRUE</CELLTEMP><CO2ABS>TRUE</CO2ABS> <CO2>TRUE</CO2><CELLPRES>TRUE</CELLPRES><IVOLT>TRUE</IVOLT><RAW>TRUE </RAW></RS232></LI840>

Тад	Parent	Value(s)	R/W	Comments
<co2></co2>	<raw></raw>	Integer	R	
<co2ref></co2ref>	<raw></raw>	Integer	R	
<h2o></h2o>	<raw></raw>	Integer	R	
<h2oref></h2oref>	<raw></raw>	Integer	R	

Tag	Parent	Value(s)	R/W	Comments
<cfg></cfg>	<li840></li840>	<outrate></outrate>		Elements within the <cfg> tag control system set-</cfg>
		<heater></heater>		tings.
		<pcomp></pcomp>		
		<doffset></doffset>		
		<filter></filter>		
		<alarms></alarms>		
		<dacs></dacs>		
		<bench></bench>		
				
		?		

<LI840><CFG><HEATER>TRUE</HEATER><PCOMP>TRUE</PCOMP></CFG></LI840>

Тад	Parent	Value(s)	R/W	Comments
<outrate></outrate>	<cfg></cfg>	0.5 to 20	R/W	Output data every N seconds (0.5 s increments).
<heater></heater>	<cfg></cfg>	TRUE FALSE	R/W	Turn heater on/off.
<pcomp></pcomp>	<cfg></cfg>	TRUE FALSE	R/W	Pressure compensation on/off.
<filter></filter>	<cfg></cfg>	0 to 20	R/W	Set a 0 to 20 second filter.
<alarms></alarms>	<cfg></cfg>	<source/> <enabled> <high> <hdead> <low> <ldead></ldead></low></hdead></high></enabled>		High and low alarm settings.
<dacs></dacs>	<cfg></cfg>	<range> <set1> <d1> <d1_0> <d1_f> <d2> <d2_0> <d2 f=""></d2></d2_0></d2></d1_f></d1_0></d1></set1></range>		DAC outputs.
<bench></bench>	<cfg></cfg>	14	R	Optical bench length (cm).

Тад	Parent	Value(s)	R/W	Comments
<enabled></enabled>	<alarms></alarms>	TRUE FALSE	R/W	Enable alarms.
<source/>	<alarms></alarms>	CO2 H2O	R/W	Channel (CO ₂ :ppm, H ₂ O:mmol/mol)
<high></high>	<alarms></alarms>	Float	R/W	High alarm on at this value.
<hdead></hdead>	<alarms></alarms>	Float	R/W	High alarm off at this value.
<low></low>	<alarms></alarms>	Float	R/W	Low alarm on at this value.
<ldead></ldead>		Float	R/W	Low alarm off at this value.

<LI840><CFG><ALARMS><ENABLED>TRUE</ENABLED><SOURCE>H2O</SOURCE> <HIGH>1600</HIGH><HDEAD>1500</HDEAD><LOW>200</LOW><LDEAD>300</LDEAD> </ALARMS></CFG></LI840>

Тад	Parent	Value(s)	R/W	Comments
<range></range>	<dacs></dacs>	2.5 5.0	R/W	Output voltage. 2.5V or 5.0V.
<set1></set1>	<dacs></dacs>	Float	R/W	Set DAC output to value (DACS must be NONE)
<set2></set2>	<dacs></dacs>	Float	R/W	Set DAC output to value (DACS must be NONE)
<d1></d1>	<dacs></dacs>	CO2 H2O H2ODEWPOINT CELLPRES CELLTEMP NONE	R/W	DAC1
<d1_0></d1_0>	<dacs></dacs>	Float	R/W	Value where DAC1 outputs 0V
<d1_f></d1_f>	<dacs></dacs>	Float	R/W	Value where DAC1 outputs full scale
<d2></d2>	<dacs></dacs>	CO2 H2O H2ODEWPOINT CELLPRES CELLTEMP NONE	R/W	DAC2
<d2_0></d2_0>	<dacs></dacs>	Float	R/W	Value where DAC2 outputs 0V
<d2_f></d2_f>	<dacs></dacs>	Float	R/W	Value where DAC2 outputs full scale

<L1840><CFG><DACS><RANGE>2.5</RANGE><D1>CO2</D1><D2>CELLTEMP</D2></DACS></CFG></L1840>

Appendix E

Tag	Parent	Value(s)	R/W	Comments
<cal></cal>	<li840></li840>	<date></date>		Calibration parameters. When calibrating,
		<h2ozero></h2ozero>		<date> must be paired with a <co2zero></co2zero></date>
		<h2ospan></h2ospan>		or a <co2span>, or a <h2ozero> or</h2ozero></co2span>
		<h2ospan2></h2ospan2>		<h2ospan>.</h2ospan>
		<h2olastzero></h2olastzero>		
		<h2olastspan></h2olastspan>		
		<h2olastspan2></h2olastspan2>		
		<h2okzero></h2okzero>		
		<h2okspan></h2okspan>		
		<h2okspan2></h2okspan2>		
		<co2zero></co2zero>		
		<co2span></co2span>		
		<co2span2></co2span2>		
		<co2lastzero></co2lastzero>		
		<co2lastspan></co2lastspan>		
		<co2lastspan2></co2lastspan2>		
		<co2kzero></co2kzero>		
		<co2kspan></co2kspan>		
		<co2kspan2></co2kspan2>		
		<poly></poly>		

Тад	Parent	Value(s)	R/W	Comments
<date></date>	<cal></cal>	20 character date	W	The date the calibration is taking place.
<h2ozero></h2ozero>	<cal></cal>	TRUE FALSE	W	Start H ₂ O zero.
<h2ospan></h2ospan>	<cal></cal>	Float (H ₂ O °C)	W	Start the first H ₂ O span.
<h2ospan2></h2ospan2>	<cal></cal>	Float (H ₂ O °C)	W	Start the second H ₂ O span.
<h2olastzero></h2olastzero>	<cal></cal>	20 character date	R	Date the LI-840A was last zeroed.
<h2olastspan></h2olastspan>	<cal></cal>	20 character date	R	Date the LI-840A was last spanned.
<h2okzero></h2okzero>	<cal></cal>	Float	R/W	Calibration constant.
<h2okspan></h2okspan>	<cal></cal>	Float	R/W	Calibration constant.
<h2okspan2></h2okspan2>	<cal></cal>	Float	R/W	Calibration constant.
<co2zero></co2zero>	<cal></cal>	TRUE FALSE	W	Start a CO_2 zero.
<co2span></co2span>	<cal></cal>	Float (CO ₂ ppm)	W	Start the first CO ₂ span.
<co2span2></co2span2>	<cal></cal>	Float (CO ₂ ppm)	W	Start the second CO ₂ span.
<co2lastzero></co2lastzero>	<cal></cal>	20 character date	R	The date the LI-840A was last zeroed.
<co2lastspan></co2lastspan>	<cal></cal>	20 character date	R	The date the LI-840A was last spanned.
<co2kzero></co2kzero>	<cal></cal>	Float	R/W	Calibration constant.
<co2kspan></co2kspan>	<cal></cal>	Float	R/W	Calibration constant.
<co2kspan></co2kspan>	<cal></cal>	Float	R/W	Calibration constant.

Appendix E

Тад	Parent	Value(s)	R/W	Comments
<poly></poly>	<li840></li840>	<li840></li840>	R/W	Serial number and factory calibration date: HGA_XXXX Day Month Year
		<bb></bb>	R/W	Band broadening.
		<xs></xs>	R/W	Cross sensitivity.
		<co2></co2>	R/W	CO ₂ coefficients.
		<h2o></h2o>	R/W	H ₂ O coefficients.
		<press></press>	R/W	Pressure correction coefficients

Examples:

To Zero:

<LI840><CAL><DATE>2015-02-07</DATE>CO2ZERO>TRUE</CO2ZERO></CAL></LI840>

To Span:

<LI840><CAL><DATE>2015-02-07</DATE><C02SPAN>1000</C02SPAN></CAL></LI840>

Possible LI-840 Replies:

<LI840><ACK>TRUE</ACK></LI840> <LI840><ERROR>{ErrorText}</ERROR></LI840>

<L1840><CAL><CO2LASTSPAN>{isodate}</CO2LASTSPAN><CO2LASTZERO>{isod-

ate}</CO2LASTZERO><CO2KZERO>{float}

</CO2KZERO><CO2KSPAN>{float}</CO2KSPAN></CAL></LI840>

Tag	Parent	Value(s)	R/W	Comments
<li840></li840>	<poly></poly>	20 character date	R/W	Serial number and factory cal- ibration date:HGA XXXX Day Month Year
<bb></bb>	<poly></poly>	Float	R/W	Band broadening value.
<xs></xs>	<poly></poly>	Float	R/W	Cross sensitivity value.
<co2></co2>	<poly></poly>	<a1><a2><a3><a4></a4></a3></a2></a1>	R/W	Floating point coefficients for the CO ₂ calibration inverse rectangular hyperbola.
<h2o></h2o>	<poly></poly>	<a1><a2><a3></a3></a2></a1>	R/W	Floating point coefficients for the H_2O calibration polynomial.
<press></press>	<poly></poly>	<a0></a0>	R/W <a1></a1>	

Тад	Parent	Value(s)	R/W	Comments
<ack></ack>	<li840></li840>	TRUE FALSE	R	Acknowledgement to a send command.
<ver></ver>	<li840></li840>	Text	R	Embedded software version.
<error></error>	<li840></li840>	Text	R	Error. Includes a message.

Warranty

Each LI-COR, inc. instrument is warranted by LI-COR, inc. to be free from defects in material and workmanship; however, LI-COR, inc.'s sole obligation under this warranty shall be to repair or replace any part of the instrument which LI-COR, inc.'s examination discloses to have been defective in material or workmanship without charge and only under the following conditions, which are:

1. The defects are called to the attention of LI-COR, inc. in Lincoln, Nebraska, in writing within one year after the shipping date of the instrument.

2. The instrument has not been maintained, repaired, or altered by anyone who was not approved by LI-COR, inc.

3. The instrument was used in the normal, proper, and ordinary manner and has not been abused, altered, misused, neglected, involved in and accident or damaged by act of God or other casualty.

4. The purchaser, whether it is a DISTRIBUTOR or direct customer of LI-COR or a DISTRIBUTOR'S customer, packs and ships or delivers the instrument to LI-COR, inc. at LI-COR inc.'s factory in Lincoln, Nebraska, U.S.A. within 30 days after LI-COR, inc. has received written notice of the defect. Unless other arrangements have been made in writing, transportation to LI-COR, inc. (by air unless otherwise authorized by LI-COR, inc.) is at customer expense.

5. No-charge repair parts may be sent at LI-COR, inc.'s sole discretion to the purchaser for installation by purchaser.

6. LI-COR, inc.'s liability is limited to repair or replace any part of the instrument without charge if LI-COR, inc.'s examination disclosed that part to have been defective in material or workmanship.

There are no warranties, express or implied, including but not limited to any implied warranty of merchantability of fitness for a particular purpose on underwater cables or on expendables such as batteries, lamps, thermocouples and calibrations.

Other than the obligation of LI-COR, inc. expressly set forth herein, LI-COR, inc. disclaims all warranties of merchantability or fitness for a particular purpose. The foregoing constitutes LI-COR, inc.'s sole obligation and liability with respect to damages resulting from the use or performance of the instrument and in no event shall LI-COR, inc. or its representatives be liable for damages beyond the price paid for the instrument, or for direct, incidental or consequential damages.

The laws of some locations may not allow the exclusion or limitation on implied warranties or on incidental or consequential damages, so the limitations herein may not apply directly. This warranty gives you specific legal rights, and you may already have other rights which vary from state to state. All warranties that apply, whether included by this contract or by law, are limited to the time period of this warranty which is a twelve-month period commencing from the date the instrument is shipped to a user who is a customer or eighteen months from the date of shipment to LI-COR, inc.'s authorized distributor, whichever is earlier.

This warranty supersedes all warranties for products purchased prior to June 1, 1984, unless this warranty is later superseded.

DISTRIBUTOR or the DISTRIBUTOR'S customers may ship the instruments directly to LI-COR if they are unable to repair the instrument themselves even though the DISTRIBUTOR has been approved for making such repairs and has agreed with the customer to make such repairs as covered by this limited warranty.

Further information concerning this warranty may be obtained by writing or telephoning Warranty manager at LI-COR, inc.

Measuring Change in a Changing World[®]



Biosciences

LI-COR, Inc. • Environmental • 4647 Superior Street • P.O. Box 4425 • Lincoln, Nebraska 68504 USA Phone: 402-467-3576 • FAX: 402-467-2819 Toll-free 1-800-447-3576 (U.S. & Canada) envsales@licor.com www.licor.com